

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Peanut R6202	Range for $z \leq 2$ according to R-Biopharm Ridascreen Fast Peanut R6202	Result (mg/kg)	Recovery (%)
FAPAS - Peanut in Garlic Powder - 27325A - May - July 2022	17.0	8.52 – 25.5	11.41	67.1
FAPAS - Peanut in Garlic Powder - 27325B - May - July 2022	NOT DETECTED	-	<1.0	-
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27319 - March - May 2022	21.0	10.5 – 31.5	29.71	141.5
FAPAS - Peanut in Chocolate - 27315A - February - March 2022	NOT DETECTED	-	<1.0	-
FAPAS - Peanut in Chocolate - 27315B - February - March 2022	20.2	10.1 – 30.3	13.37	66.2
FAPAS - Peanut in Garlic Powder - 27297A - June - August 2021	97.0	48.5 – 146	72.4	74.6
FAPAS - Peanut in Garlic Powder	NOT DETECTED	-	<1.0	-

27297B - June - August 2021				
-----------------------------	--	--	--	--

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS - Allergens on Swab - 27383A - January - March 2024	DETECTED	DETECTED	69.0
FAPAS - Allergens on Swab - 27383A - January - March 2024	NOT DETECTED	NOT DETECTED	94.0
FAPAS - Nuts In Chocolate - 27312 - November 2021 - February 2022	DETECTED	DETECTED	98.0

Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27353 - March - May 2023	15.3	15.0	-0.1