


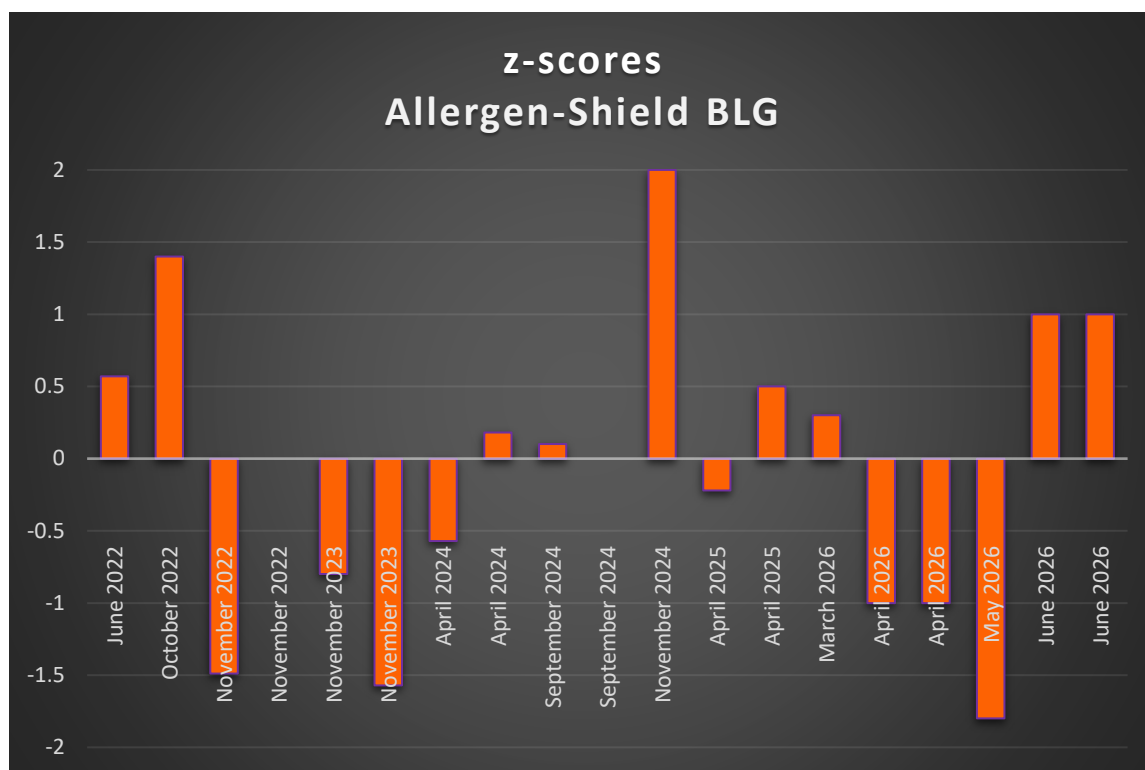

Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1077 – June 2026	0.3	0.4	1.00
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1177 – June 2026	NOT DETECTED	<0.15	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1277 – June 2026	0.5	0.7	1.00
FAPAS Milk & Gluten in Infant Soya Formula 27463A March – May 2026	NOT DETECTED	<0.15	-
FAPAS Milk & Gluten in Infant Soya Formula 27463B March – May 2026	2.72	1.50	-1.8

BIPEA Food Allergens - Milk 77C - Infant Formula – 23 - 0777 – April 2026	0.4	0.2	-1.0
BIPEA Food Allergens - Milk 77C - Infant Formula – 23 - 0877 – April 2026	NOT DETECTED	<0.15	-
BIPEA Food Allergens - Milk 77C - Infant Formula – 22 - 0977 – April 2026	0.9	0.6	-1.0
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27459 January – March 2026	2.06	2.22	0.3
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 21 - 0777 – April 2025	NOT DETECTED	<0.1	-
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 21 - 0877 – April 2025	1.3	1.2	-0.22
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 20 - 0977 – April 2025	0.5	0.6	0.50
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 20 - 0777 – November 2024	NOT DETECTED	<0.1	-
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 20 - 0877 – November 2024	NO VALUE WAS ASSIGNED	<0.1	-
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 19 - 0977 – November 2024	0.4	0.7	2.0
FAPAS Gluten, Egg and Milk in Cake Mix 27400A July – September 2024	2.91	2.97	0.1
FAPAS Gluten, Egg and Milk in Cake Mix 27400B July – September 2024	7.57	7.63	0.0
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula – 19 - 0777 – April 2024	1.4	1.2	-0.57
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula – 19 - 0877 – April 2024	2.1	2.2	0.18

BIPEA Food Allergens - Milk 77C - 18 - Infant Formula – 18 - 0977 – April 2024	NOT DETECTED	<0.1	-
BIPEA Allergens in Wines 71 - 32 - White Wine – 05 - 1571 – February 2024	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1077 – January 2024	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1177 – January 2024	NOT DETECTED	<0.1	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1277 – January 2024	DETECTED	DETECTED	-
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula – 18 - 0777 – November 2023	2.4	1.3	-1.57
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula – 18 - 0877 – November 2023	0.6	0.4	-0.8
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula – 17 - 0977 – November 2023	NOT DETECTED	<0.1	-
BIPEA Food Allergens - Milk 77C - 15 - 0204 Infant Formula – 16 - 0777 – November 2022	0.6	0.6	0.00
BIPEA Food Allergens - Milk 77C - 15 - 0204 Infant Formula – 17 - 0877 – November 2022	6.1	2.9	-1.49
BIPEA Food Allergens - Milk 77C - 15 - 0204 Infant Formula – 15 - 0977 – November 2022	NOT DETECTED	<0.1	-
BIPEA Allergens in Wines 71 - 28 - Rose Wine - 04 - 1471 – October 2022	0.28	0.42	1.40
BIPEA Allergens in Industrial Process Water 77D – 0894 Industrial Process Water - 13 – 1077 – June 2022	NOT DETECTED	<0.1	-
BIPEA Allergens in Industrial Process Water 77D - 0894 Industrial Process Water - 13 – 1177 – June 2022	DETECTED	DETECTED	-

BIPEA Allergens in Industrial Process Water 77D - 13 2021-2022 - 0894 Industrial Process Water - 13 - 1277 - June 2022	DETECTED	DETECTED	-
BIPEA Allergens in Wines 71 - 27 - 0902 - Red Wine - 03 - 1671 - June 2022	0.21	0.25	0.57

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Milk in Infant Breakfast Cereal 27435A June - August 2025	DETECTED	DETECTED	-
FAPAS Milk in Infant Breakfast Cereal 27435B June - August 2025	NOT DETECTED	<0.15	-
FAPAS Milk & Gluten in Infant Soya Formula 27422A February - March 2025	DETECTED	DETECTED	-
FAPAS Milk & Gluten in Infant Soya Formula 27422B February - March 2025	NOT DETECTED	<0.15	-
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27418 January - March 2025	DETECTED	DETECTED	-
FAPAS Milk in Infant Breakfast Cereal - 27398A - June - August 2024	NOT DETECTED	<0.15	-
FAPAS Milk in Infant Breakfast Cereal - 27398B - June - August 2024	DETECTED	DETECTED	-
FAPAS Milk and Gluten in Infant Soya Formula 27386A February - April 2024	NOT DETECTED	<0.15	-
FAPAS Milk and Gluten in Infant Soya Formula 27386B February - April 2024	DETECTED	DETECTED	-

FAPAS Allergens in Cooked Biscuit 27382 January – March 2024	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 19 - 1077 – January 2025	DETECTED	DETECTED	100.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 18 - 1177 – January 2025	NOT DETECTED	<0.15	100.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 18 - 1277 – January 2025	DETECTED	DETECTED	100.0

Proficiency Test	Assigned value (mg/kg) according to R- Biopharm Ridascreen Fast β-Lactoglobulin R4912	Range for $z \leq 2$ according to R- Biopharm Ridascreen Fast Fast β- Lactoglobulin R4912	Result (mg/kg)	Recovery (%)
FAPAS Gluten, Milk & Egg in Cake Mix 27438A July – August 2025	2.12	1.06 – 3.18	2.55	120.3
FAPAS Gluten, Milk & Egg in Cake Mix 27438B July – August 2025	3.44	1.72 – 5.16	5.14	149.4
FAPAS Milk & Gluten in Plant Based Milk Alternative 27445A September – October 2025	NOT DETECTED	-	<0.15	-
FAPAS Milk & Gluten in Plant Based Milk Alternative 27445B September – October 2025	3.20	1.60 – 4.80	2.80	87.5
FAPAS Allergens in Cake Mix - 27364A - July – September 2023	2.30	1.15 – 3.45	2.21	96.1

FAPAS Allergens in Cake Mix - 27364B - July – September 2023	3.70	1.85 – 5.55	3.90	105.4
FAPAS Milk in Infant Breakfast Cereal - 27362A - July – August 2023	NOT DETECTED	-	<0.15	-
FAPAS Milk in Infant Breakfast Cereal - 27362B - July – August 2023	1.60	0.800 – 2.40	0.96	60.0
FAPAS Allergens in Cake Mix - 27331A - July – September 2022	1.00	0.500 – 1.50	1.31	131.0
FAPAS Allergens in Cake Mix - 27331B - July – September 2022	0.600	0.300 – 0.900	0.656	109.3
FAPAS Milk in Infant Breakfast Cereal - 27329A - June – August 2022	1.11	0.555 – 1.67	1.63	146.8
FAPAS Milk in Infant Breakfast Cereal - 27329B - June – August 2022	NOT DETECTED	-	<0.15	-
FAPAS Gluten & Milk in Infant Soya Formula - 27316A - February – April 2022	DETECTED	-	DETECTED	-
FAPAS Gluten & Milk in Infant Soya Formula - 27316B - February – April 2022	NOT DETECTED	-	<0.15	-
FAPAS Allergens in Cooked Biscuit – 27314 - January – March 2022	1.51	0.752 – 2.26	1.51	100.0
FAPAS Milk in Infant Soya Formula - 27313A - January – February 2022	NOT DETECTED	-	<0.15	-
FAPAS Milk in Infant Soya Formula - 27313B - January – February 2022	DETECTED	-	DETECTED	-
FAPAS Milk in Infant Breakfast Cereal - 27322A - September – October 2021	NOT DETECTED	-	<0.15	-
FAPAS Milk in Infant Breakfast Cereal - 27322B - September – October 2021	1.52	0.76 – 2.28	1.35	88.8



FAPAS Milk in Infant Formula (hydrolyzed) - 27304A - August - October 2021	NOT DETECTED	-	<0.15	-
FAPAS Milk in Infant Formula (hydrolyzed) - 27304B - August - October 2021	DETECTED	-	DETECTED	-
FAPAS Allergens in Cake Mix - 27301A - July - September 2021	1.19	0.595 - 1.790	0.89	74.8
FAPAS Allergens in Cake Mix - 27301B - July - September 2021	0.545	0.273 - 0.818	0.47	86.2
FAPAS Allergens in Cooked Biscuit - 27286 - January - March 2021	3.58	1.79 - 5.37	3.06	85.5