


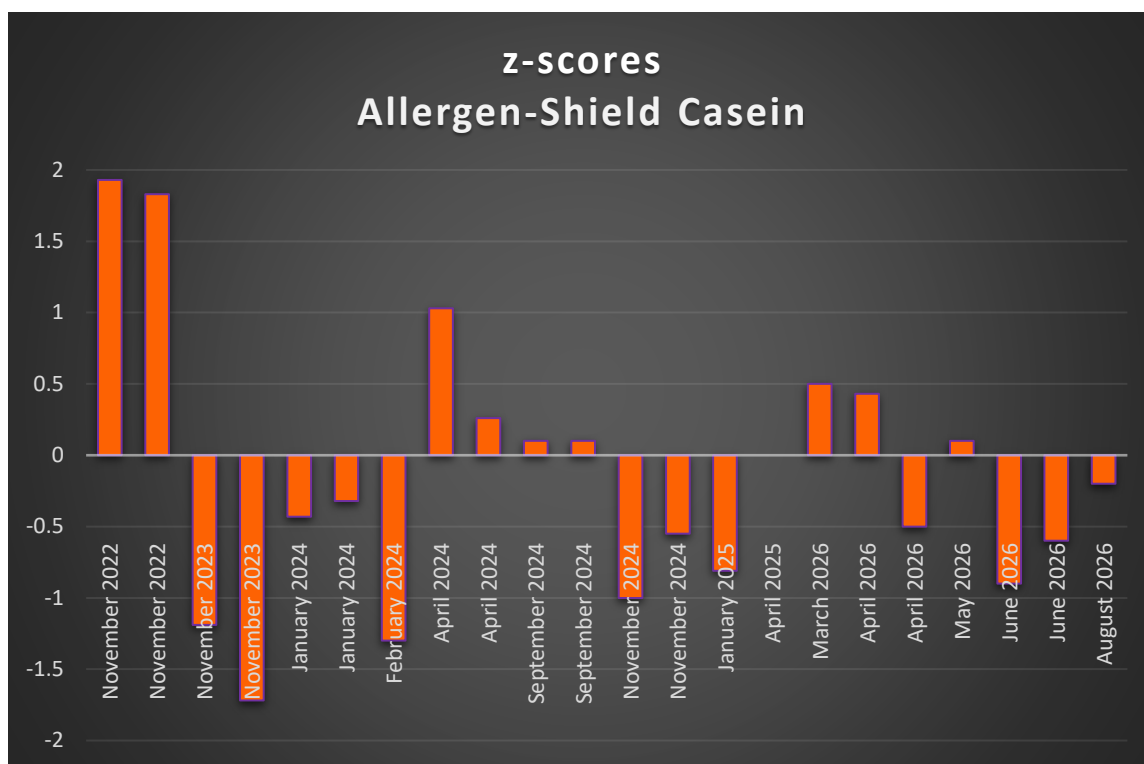

Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS Milk in Infant Breakfast Cereal 27475A June – August 2026	18.1	17.1	-0.2
FAPAS Milk in Infant Breakfast Cereal 27475B June – August 2026	NOT DETECTED	<0.5	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1077 – June 2026	17.8	11.6	-0.90
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1177 – June 2026	NOT DETECTED	<0.5	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1277 – June 2026	24.8	20.4	-0.60
FAPAS Milk & Gluten in Infant Soya Formula 27463A March – May 2026	NOT DETECTED	<0.5	-
FAPAS Milk & Gluten in Infant Soya Formula 27463B March – May 2026	18.1	18.7	0.1

BIPEA Food Allergens - Milk 77C - Infant Formula – 23 - 0777 – April 2026	6.7	7.7	0.43
BIPEA Food Allergens - Milk 77C - Infant Formula – 23 - 0877 – April 2026	NOT DETECTED	<0.5	-
BIPEA Food Allergens - Milk 77C - Infant Formula – 22 - 0977 – April 2026	20.2	16.7	-0.5
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27459 January – March 2026	19.9	22.5	0.5
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 21 - 0777 – April 2025	NOT DETECTED	<0.5	-
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 21 - 0877 – April 2025	26.6	DETECTED	-
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 20 - 0977 – April 2025	10.1	10.1	0.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 19 - 1077 – January 2025	9.9	7.8	-0.81
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 18 - 1177 – January 2025	NOT DETECTED	<0.5	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 18 - 1277 – January 2025	DETECTED	DETECTED	-
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 20 - 0777 – November 2024	NOT DETECTED	<0.5	-
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 20 - 0877 – November 2024	2.0	1.6	-1.0
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 19 - 0977 – November 2024	15.7	13.4	-0.55
FAPAS Gluten, Egg and Milk in Cake Mix 27400A July – September 2024	14.4	14.7	0.1
FAPAS Gluten, Egg and Milk in Cake Mix 27400B July – September 2024	30.4	31.2	0.1

BIPEA Food Allergens - Milk 77C - 18 - Infant Formula – 19 - 0777 – April 2024	20.5	25.8	1.03
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula – 19 - 0877 – April 2024	33.6	35.8	0.26
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula – 18 - 0977 – April 2024	NOT DETECTED	<0.5	-
BIPEA Allergens in Wines 71 - 32 - White Wine – 11 - 0671 – February 2024	0.51	0.36	-1.30
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1077 – January 2024	30.8	26.9	-0.32
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1177 – January 2024	NOT DETECTED	<1.0	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1277 – January 2024	34.0	27.6	-0.43
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula – 18 - 0777 – November 2023	8.0	4.3	-1.72
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula – 18 - 0877 – November 2023	4.1	2.2	-1.19
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula – 17 - 0977 – November 2023	NOT DETECTED	<1.0	-
BIPEA Food Allergens - Milk 77C - 15 - 0204 Infant Formula – 16 - 0777 – November 2022	5.4	8.6	1.83
BIPEA Food Allergens - Milk 77C - 15 - 0204 Infant Formula – 17 - 0877 – November 2022	27.6	44.9	1.93
BIPEA Food Allergens - Milk 77C - 15 - 0204 Infant Formula – 15 - 0977 – November 2022	NOT DETECTED	<2.5	-
BIPEA Allergens in Wines 71 - 28 - Rose Wine – 09 - 1271 – October 2022	0.30	<0.60	-
BIPEA Allergens in Industrial Process Water 77D – 0894 Industrial Process Water - 13 – 1077 – June 2022	NOT DETECTED	<2.5	-

BIPEA Allergens in Industrial Process Water 77D - 0894 Industrial Process Water - 13 - 1177 - June 2022	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - 13 2021- 2022 - 0894 Industrial Process Water - 13 -1277 - June 2022	DETECTED	DETECTED	-
BIPEA Allergens in Wines 71 - 27 - 0902 - Red Wine - 08 - 0971- June 2022	NOT DETECTED	<0.25	-

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Casein R4612	Range for $z \leq 2$ according to R- Biopharm Ridascreen Fast Casein R4612	Result (mg/kg)	Recovery (%)
FAPAS Gluten, Milk & Egg in Cake Mix 27438A July - August 2025	17.5	8.74 - 26.2	20.8	118.9
FAPAS Gluten, Milk & Egg in Cake Mix 27438B July - August 2025	26.5	13.3 - 39.8	39.6	149.4
FAPAS Milk in Infant Breakfast Cereal 27435A June - August 2025	14.0	7.01 - 21.0	13.7	97.9
FAPAS Milk in Infant Breakfast Cereal 27435B June - August 2025	NOT DETECTED	-	<0.50	-
FAPAS Milk & Gluten in Infant Soya Formula 27422A February - March 2025	19.7	9.86 - 29.6	22.2	112.7
FAPAS Milk & Gluten in Infant Soya Formula 27422B February - March 2025	NOT DETECTED	-	<0.50	-
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27418 January - March 2025	17.9	8.93 - 26.8	17.8	99.4

FAPAS Milk in Infant Breakfast Cereal - 27398A - June – August 2024	NOT DETECTED	-	<0.5	-
FAPAS Milk in Infant Breakfast Cereal - 27398B - June – August 2024	19.1	9.55 – 28.6	13.4	70.2
FAPAS Milk and Gluten in Infant Soya Formula 27386A February – April 2024	NOT DETECTED	-	<0.5	-
FAPAS Milk and Gluten in Infant Soya Formula 27386B February – April 2024	13.3	6.65 – 20.0	17.7	133.1
FAPAS Allergens in Cooked Biscuit 27382 January – March 2024	13.0	6.50 – 19.5	15.02	115.5
FAPAS Allergens in Cake Mix - 27364A - July – September 2023	14.2	7.09 – 21.3	14.1	99.3
FAPAS Allergens in Cake Mix - 27364B - July – September 2023	35.0	17.5 – 52.5	31.4	89.7
FAPAS Milk in Infant Breakfast Cereal - 27362A - July – August 2023	NOT DETECTED	-	<1.0	-
FAPAS Milk in Infant Breakfast Cereal - 27362B - July – August 2023	18.0	9.00 – 27.0	16.1	89.4
FAPAS Allergens in Cake Mix - 27331A - July – September 2022	12.4	6.18 – 18.5	10.75	86.7
FAPAS Allergens in Cake Mix - 27331B - July – September 2022	5.63	2.82 – 8.45	4.76	84.5
FAPAS Milk in Infant Breakfast Cereal - 27329A - June – August 2022	15.2	7.58 – 22.7	20.1	132.2

FAPAS Milk in Infant Breakfast Cereal - 27329B - June – August 2022	NOT DETECTED	-	<1.0	-
FAPAS Gluten & Milk in Infant Soya Formula - 27316A - February – April 2022	31.0	15.5 – 46.5	23.04	74.3
FAPAS Gluten & Milk in Infant Soya Formula - 27316B - February – April 2022	NOT DETECTED	-	<1.0	-
FAPAS Allergens in Cooked Biscuit - 27314 - January – March 2022	DETECTED	-	DETECTED	-
FAPAS Milk in Infant Breakfast Cereal - 27322A - September – October 2021	NOT DETECTED	-	<1.0	-
FAPAS Milk in Infant Breakfast Cereal - 27322B - September – October 2021	21.1	10.5 – 31.6	19.59	92.8
FAPAS Milk in Infant Formula (hydrolyzed) - 27304A - August – October 2021	NOT DETECTED	-	<1.0	-
FAPAS Milk in Infant Formula (hydrolyzed) - 27304B - August – October 2021	DETECTED	-	DETECTED	-
FAPAS Allergens in Cake Mix - 27301A - July – September 2021	11.0	5.5 – 16.5	13.99	127.2
FAPAS Allergens in Cake Mix - 27301B - July – September 2021	6.0	3.0 – 9.0	5.2	86.7
FAPAS Allergens in Cooked Biscuit - 27286 - January – March 2021	30.1	15.0 – 45.1	28.7	95.3

