


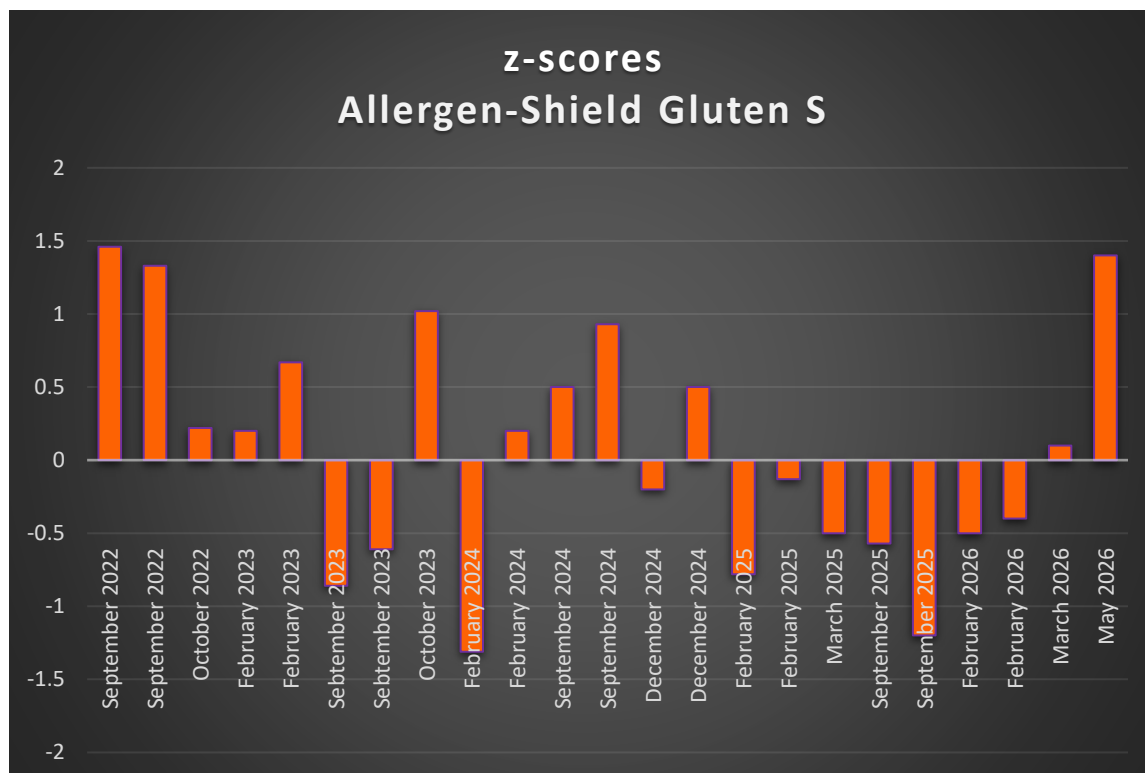

Proficiency Test	Assigned value (mg/L or mg/kg)	Result (mg/L or mg/kg)	Z-score
FAPAS Milk & Gluten in Infant Soya Formula 27463A March – May 2026	NOT DETECTED	<5.0	-
FAPAS Milk & Gluten in Infant Soya Formula 27463B March – May 2026	17.5	23.8	1.4
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27459 January – March 2026	11.3	11.7	0.1
BIPEA Food Allergens 77A - Gluten – Cake 8 - 3477 – February 2026	46.0	39.0	-0.5
BIPEA Food Allergens 77A - Gluten – Cake 8 - 3577 – February 2026	17.0	15.0	-0.4

BIPEA Food Allergens 77A - Gluten – Cake 8 - 3677 – February 2026	NOT DETECTED	<5.0	-
BIPEA Food Allergens 77A - Gluten – Infant Flour 14 - 0177 – September 2025	73.0	60.0	-0.57
BIPEA Food Allergens 77A - Gluten – Infant Flour 12 - 0277 – September 2025	NOT DETECTED	<5.0	-
BIPEA Food Allergens 77A - Gluten – Infant Flour 13 - 0377 – September 2025	11.0	5.00	-1.2
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27418 January – March 2025	22.2	19.3	-0.5
BIPEA Food Allergens 77A - Gluten – Cake 7 - 3477 – February 2025	NOT DETECTED	<5.0	-
BIPEA Food Allergens 77A - Gluten – Cake 7 - 3577 – February 2025	32.0	25.0	-0.78
BIPEA Food Allergens 77A - Gluten – Cake 7 - 3677 – February 2025	66.0	64.0	-0.13
FAPAS Gluten, Egg and Milk in Cake Mix 27400A July – September 2024	12.8	12.2	-0.2
FAPAS Gluten, Egg and Milk in Cake Mix 27400B July – September 2024	24.3	27.4	0.5
BIPEA Food Allergens 77A - Gluten – Infant Flour 13 - 0177 - September 2024	24.0	28.0	0.5
BIPEA Food Allergens 77A - Gluten – Infant Flour 11 - 0277 - September 2024	79.0	92.0	0.93
BIPEA Food Allergens 77A - Gluten – Infant Flour 12 - 0377 - September 2024	NOT DETECTED	<5.0	-
BIPEA Food Allergens 77A - Gluten - Cake 06 - 3477 - February 2024	24.0	25.0	0.20

BIPEA Food Allergens 77A - Gluten - Cake 06 - 3577 - February 2024	NOT DETECTED	<5.0	-
BIPEA Food Allergens 77A - Gluten - Cake 06 - 3677 - February 2024	54.0	37.0	-1.31
BIPEA Allergens in Wines 71 - Rose Wine – 04 - 1771 – October 2023	11.3	15.1	1.02
BIPEA Food Allergens Gluten 77A - Infant Flour – 12 - 0177 - September 2023	NOT DETECTED	<5.0	-
BIPEA Food Allergens Gluten 77A - Infant Flour – 10 - 0277 - September 2023	53.0	39.0	-0.61
BIPEA Food Allergens Gluten 77A - Infant Flour – 11 - 0377 - September 2023	16.0	10.0	-0.86
BIPEA Food Allergens Gluten 77A - Cake – 05 - 3477 - February 2023	66.0	72.0	0.67
BIPEA Food Allergens Gluten 77A - Cake – 05 - 3577 - February 2023	33.0	34.0	0.2
BIPEA Food Allergens Gluten 77A - Cake – 05 - 3677 – February 2023	NOT DETECTED	<5.0	-
BIPEA Allergens in Wines 71 - Rose Wine – 03 - 1771 – October 2022	10.33	11.22	0.22
BIPEA Food Allergens Gluten 77A - Infant Flour – 11 - 0177 - September 2022	21.0	33.0	1.33
BIPEA Food Allergens Gluten 77A - Infant Flour - 9 - 0277- September 2022	NOT DETECTED	<5.0	-
BIPEA Food Allergens Gluten 77A - Infant Flour – 10 - 0377 - September 2022	72.0	91.0	1.46
BIPEA Allergens in Wines 71 - Red Wine – 02 - 1971 – June 2022	NOT DETECTED	<10.0	-

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Allergens on Cotton Swabs 27440A August – October 2025	DETECTED	DETECTED	100.0
FAPAS Allergens on Cotton Swabs 27440B August – October 2025	NOT DETECTED	NOT DETECTED	100.0
FAPAS Gluten, Egg and Milk on Swabs (Cotton) 27366A September – November 2023	NOT DETECTED	NOT DETECTED	91.0
FAPAS Gluten, Egg and Milk on Swabs (Cotton) 27366B September – November 2023	DETECTED	DETECTED	97.0

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Gliadin R7001	Range for $ z \leq 2$ according to R-Biopharm Ridascreen Gliadin R7001	Result (mg/kg)	Recovery (%)
FAPAS Gluten & Sesame in Spice - Cumin 27464A February – April 2026	31.4	15.7 – 47.2	25.2	80.3
FAPAS Gluten & Sesame in Spice - Cumin 27464B February – April 2026	NOT DETECTED	-	<5.0	-
FAPAS Gluten, Milk & Egg in Cake Mix 27438A July – August 2025	20.5	10.3 – 30.8	12.9	62.9
FAPAS Gluten, Milk & Egg in Cake Mix 27438B July – August 2025	38.9	19.4 – 58.3	35.0	90.0
FAPAS Milk & Gluten in Plant Based Milk Alternative 27445A September – October 2025	NOT DETECTED	-	<5.0	-

FAPAS Milk & Gluten in Plant Based Milk Alternative 27445B September – October 2025	23.0	11.5 – 34.5	25.0	108.7
FAPAS Gluten in Oat Based Foodstuff 27439A July – September 2025	NOT DETECTED	-	<5.0	-
FAPAS Gluten in Oat Based Foodstuff 27439B July – September 2025	24.4	12.2 – 36.7	20.8	85.2
FAPAS Milk & Gluten in Infant Soya Formula 27422A February – March 2025	37.5	18.7 – 56.2	25.6	68.3
FAPAS Milk & Gluten in Infant Soya Formula 27422B February – March 2025	NOT DETECTED	-	<5.0	-
FAPAS Allergens in Sausage Meat (Raw) 27427A March – May 2025	18.5	9.23 – 27.7	20.57	111.2
FAPAS Allergens in Sausage Meat (Raw) 27427B March – May 2025	NOT DETECTED	-	<5.0	-
FAPAS Gluten & Sesame in Cumin 27423A February – April 2025	50.1	25.1 – 75.2	49.5	98.8
FAPAS Gluten & Sesame in Cumin 27423B February – April 2025	NOT DETECTED	-	<5.0	-
FAPAS Milk and Gluten in Infant Soya Formula 27386A February – April 2024	NOT DETECTED	-	<5.0	-
FAPAS Milk and Gluten in Infant Soya Formula 27386B February – April 2024	23.1	11.5 – 34.6	16.8	72.7

FAPAS Allergens in Cooked Biscuit 27382 January – March 2024	38.1	19.1 – 57.2	46.0	120.7
FAPAS Allergens in Cake Mix 27377A November 2023 – January 2024	10.5	5.27 – 15.8	14.7	140.0
FAPAS Allergens in Cake Mix 27377B November 2023 – January 2024	39.3	19.7 – 59.0	43.8	111.5
FAPAS Allergens in Cake Mix 27364A July – September 2023	18.9	9.44 – 28.3	16.7	88.4
FAPAS Allergens in Cake Mix 27364B July – September 2023	41.6	20.8 – 62.4	53.0	127.4
FAPAS Allergens in Cake Mix 27331A July – September 2022	13.9	6.94 – 20.8	12.62	90.8
FAPAS Allergens in Cake Mix 27331B July – September 2022	5.53	2.76 – 8.29	5.77	104.3
FAPAS Gluten in Oat Based Foodstuff 27332A July – September 2022	NOT DETECTED	-	<5.0	-
FAPAS Gluten in Oat Based Foodstuff 27332B July – September 2022	22.7	11.4 – 34.1	26.61	117.2
FAPAS Gluten in Sausage Meat (Raw) 27321A March – May 2022	DETECTED	-	DETECTED	-
FAPAS Gluten in Sausage Meat (Raw) 27321B March – May 2022	NOT DETECTED	-	<5.0	-
FAPAS Gluten & Sesame in Cumin 27317A February – April 2022	18.7	9.36 – 28.1	22.28	119.1

FAPAS Gluten & Sesame in Cumin 27317B February – April 2022	NOT DETECTED	-	<5.0	-
FAPAS Gluten & Milk in Infant Soya Formula 27316A February – April 2022	20.8	10.4 – 31.1	27.98	134.5
FAPAS Gluten & Milk in Infant Soya Formula 27316B February – April 2022	NOT DETECTED	-	<5.0	-
FAPAS Allergens in Cooked Biscuit 27314 January – March 2022	56.3	28.1 – 84.4	60.1	106.7
FAPAS Allergens in Cake Mix 27311A November 2021 – February 2022	21.1	10.6 – 31.7	20.48	97.1
FAPAS Allergens in Cake Mix 27311B November 2021 – February 2022	10.3	5.13 – 15.4	9.35	90.8
FAPAS Allergens in Cake Mix 27301A July – September 2021	8.11	4.06 – 12.20	8.68	107.0
FAPAS Allergens in Cake Mix 27301B July – September 2021	16.0	7.99 – 24.00	24.05	150.3
FAPAS Gluten in Oat Based Foodstuff 27302A July – September 2021	16.8	8.39 – 25.20	18.63	110.9
FAPAS Gluten in Oat Based Foodstuff 27302B July – September 2021	NOT DETECTED	-	<5.0	-
FAPAS Gluten & Sesame in Cumin 27289A February – April 2021	NOT DETECTED	-	<5.0	-
FAPAS Gluten & Sesame in Cumin 27289B February – April 2021	20.3	10.1 – 30.4	21.09	103.9

FAPAS Allergens in Cooked Biscuit 27286 January – March 2021	35.6	17.8 – 53.3	35.8	100.6
FAPAS Gluten in Cake Mix 27271A May – July 2020	39.2	19.6 – 58.8	43.3	110.5
FAPAS Gluten in Cake Mix 27271B May – July 2020	NOT DETECTED	-	<5.0	-

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Allergens on Cotton Swabs 27440A August – October 2025	DETECTED	DETECTED	100.0
FAPAS Allergens on Cotton Swabs 27440A August – October 2025	NOT DETECTED	NOT DETECTED	100.0
FAPAS Gluten, Egg and Milk on Swabs (Cotton) 27366A September – November 2023	NOT DETECTED	NOT DETECTED	91.0
FAPAS Gluten, Egg and Milk on Swabs (Cotton) 27366B September – November 2023	DETECTED	DETECTED	97.0

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Gliadin Competitive R7021	Range for $ z \leq 2$ according to R-Biopharm Ridascreen Gliadin Competitive R7021	Result (mg/kg)	Recovery (%)
FAPAS Gluten in Beer 27308A October – December 2022	NOT DETECTED	-	<5.0	-
FAPAS Gluten in Beer 27308B October – December 2022	17.0	8.5 – 25.5	17.2	101.2

FAPAS Gluten in Beer 27308A October – December 2021	12.2	6.09 – 18.3	14.6	119.7
FAPAS Gluten in Beer 27308B October – December 2021	NOT DETECTED	-	<10	-