

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm RIDASCREEN FAST Hazelnut R6802	Range for $ z \leq 2$ according to R-Biopharm RIDASCREEN FAST Hazelnut R6802	Result (mg/kg)	Recovery (%)
FAPAS - Hazelnut in Chocolate - 27327A May - July 2022	NOT DETECTED	-	<0.25	-
FAPAS - Hazelnut in Chocolate - 27327B May - July 2022	10.3	5.13 - 15.4	15.3	148.5
FAPAS - Hazelnut and Peanut in Cooked Biscuit -27319 - March - May 2022	18.0	9.01 - 27.0	17.69	98.3
FAPAS - Hazelnut in Chocolate - 27299A - May - July 2021	16.4	8.19 - 24.6	11.22	68.4
FAPAS - Hazelnut in Chocolate - 27299B - May - July 2021	NOT DETECTED	-	<0.25	-
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27291 March - May 2021	10.8	5.42 - 16.3	12.3	113.9

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS - Nuts In Chocolate - 27378 - November 2023 - January 2024	NOT DETECTED	<0.25	100.0
FAPAS - Nuts In Chocolate - 27343 -November 2022 - January 2023	DETECTED	DETECTED	100.0
FAPAS - Nuts In Chocolate - 27312 -November 2021 - February 2022	NOT DETECTED	<0.25	96.0

Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27353 March - May 2023	6.76	7.16	0.2
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 - 0795 - May 2023 - Cake - 11-2877	58.7	46.8	-0.61
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 - 0795 - May 2023 - Cake - 09-2977	32.7	13.8	-1.84
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 - 0795 - May 2023 - Cake - 13-3077	NOT DETECTED	<0.25	-