



Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March - April 2026	3.69 (mg/kg of allergenic protein)	4.01	0.3
	22.1	26.8	0.9
BIPEA - Food Allergens - Nuts 77G - 19 - May 2025 - Cake - 13-2877	532.9	667.4	0.77
BIPEA - Food Allergens - Nuts 77G - 19 - May 2025 - Cake - 11-2977	318.6	417.3	0.84
BIPEA - Food Allergens - Nuts 77G - 19 - May 2025 - Cake - 15-3077	NOT DETECTED	<1.0	-

FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27353 March - May 2023	6.76	7.16	0.2
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 - 0795 - May 2023 - Cake - 11-2877	58.7	46.8	-0.61
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 - 0795 - May 2023 - Cake - 09-2977	32.7	13.8	-1.84
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 - 0795 - May 2023 - Cake - 13-3077	NOT DETECTED	<1.0	-

<b>Proficiency Test</b>	<b>Assigned value (mg/kg) according to R-Biopharm RIDASCREEN FAST Hazelnut R6802</b>	<b>Range for <math> z  \leq 2</math> according to R- Biopharm RIDASCREEN FAST Hazelnut R6802</b>	<b>Result (mg/kg)</b>	<b>Recovery (%)</b>
FAPAS - Hazelnut in Chocolate - 27327A May - July 2022	NOT DETECTED	-	<1.0	-
FAPAS - Hazelnut in Chocolate - 27327B May - July 2022	10.3	5.13 - 15.4	15.3	148.5
FAPAS - Hazelnut and Peanut in Cooked Biscuit -27319 - March - May 2022	18.0	9.01 - 27.0	17.69	98.3
FAPAS - Hazelnut in Chocolate - 27299A - May - July 2021	16.4	8.19 - 24.6	11.22	68.4
FAPAS - Hazelnut in Chocolate - 27299B - May - July 2021	NOT DETECTED	-	<1.0	-
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27291 March - May 2021	10.8	5.42 - 16.3	12.3	113.9

<b>Proficiency Test</b>	<b>Assignment</b>	<b>Result (mg/kg)</b>	<b>Satisfactory results / results agreeing with assignment (%)</b>
FAPAS - Nuts In Chocolate - 27453 - November 2025 – January 2026	NOT DETECTED	<1.0	97.0
FAPAS - Hazelnut in Chocolate-27433A - May – July 2025	DETECTED	DETECTED	100.0
FAPAS - Hazelnut in Chocolate-27433B - May – July 2025	NOT DETECTED	<1.0	95.0
FAPAS - Hazelnut and Peanut in Cooked Biscuit -27425 - March – April 2025	DETECTED	DETECTED	95.0
BIPEA - Food Allergens - Nuts 77G - 17 - May 2024 - Cake - 12-2877	DETECTED	DETECTED	100.0
BIPEA - Food Allergens - Nuts 77G - 17 - May 2024 - Cake - 10-2977	NOT DETECTED	<1.0	75.0
BIPEA - Food Allergens - Nuts 77G - 17 - May 2024 - Cake - 14-3077	DETECTED	DETECTED	100.0
FAPAS – Hazelnut and Peanut in Cooked Biscuit - 27389 - March – May 2024	DETECTED	5.70	96.0
FAPAS - Nuts In Chocolate - 27378 - November 2023 – January 2024	NOT DETECTED	<1.0	100.0
FAPAS - Nuts In Chocolate - 27343 -November 2022 – January 2023	DETECTED	DETECTED	100.0
FAPAS - Nuts In Chocolate - 27312 -November 2021 – February 2022	NOT DETECTED	<1.0	96.0