


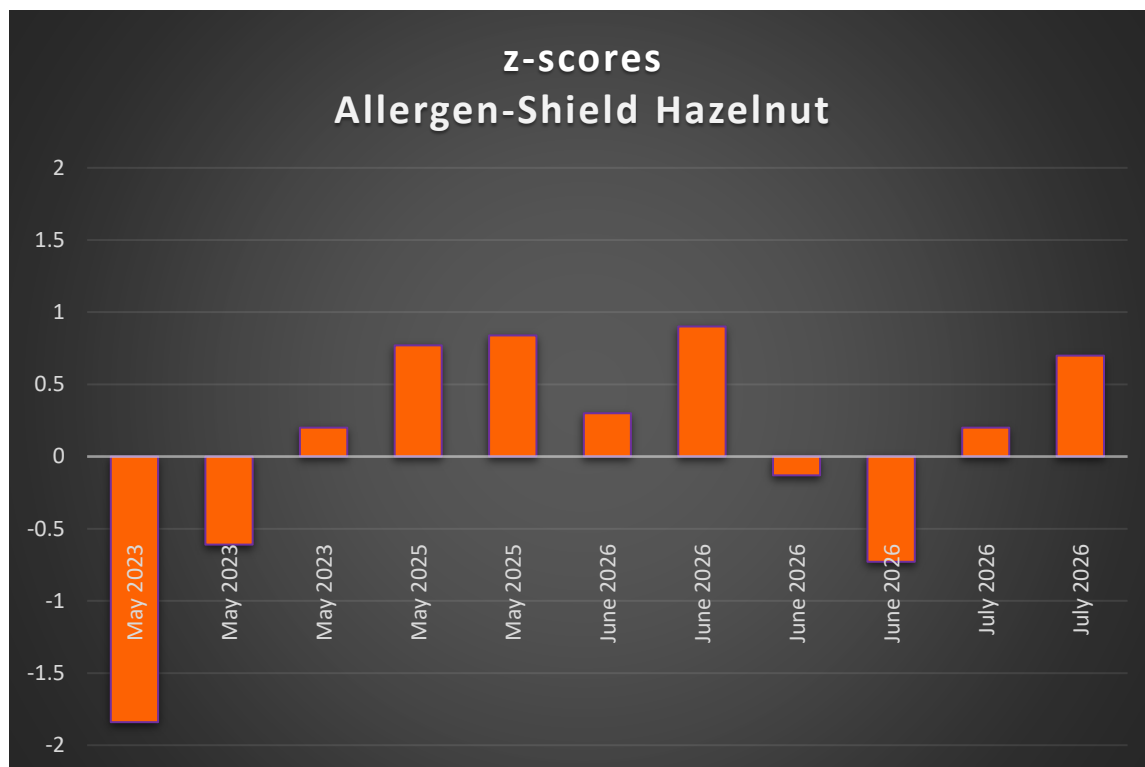

Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS - Hazelnut in Chocolate – 27473A May – July 2026	NOT DETECTED	<1.0	-
FAPAS - Hazelnut in Chocolate – 27473B May – July 2026	3.32 (mg/kg of allergenic protein)	3.90	0.7
	24.5	26.0	0.2
BIPEA - Food Allergens - Nuts 77G - 21 - May 2026 - Cake - 14-2877	NOT DETECTED	<1.0	-
BIPEA - Food Allergens - Nuts 77G - 21 - May 2026 - Cake - 12-2977	59.0	57.1	-0.13

BIPEA - Food Allergens - Nuts 77G - 21 - May 2026 - Cake - 16-3077	50.5	41.3	-0.73
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March – April 2026	3.69 (mg/kg of allergenic protein)	4.01	0.3
	22.1	26.8	0.9
BIPEA - Food Allergens - Nuts 77G – 19 - May 2025 - Cake - 13-2877	532.9	667.4	0.77
BIPEA - Food Allergens - Nuts 77G – 19 - May 2025 - Cake - 11-2977	318.6	417.3	0.84
BIPEA - Food Allergens - Nuts 77G – 19 - May 2025 - Cake - 15-3077	NOT DETECTED	<1.0	-
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27353 March – May 2023	6.76	7.16	0.2
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 – 0795 - May 2023 - Cake - 11-2877	58.7	46.8	-0.61
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 – 0795 - May 2023 - Cake - 09-2977	32.7	13.8	-1.84
BIPEA - Food Allergens - Nuts 77G - 15 2022-2023 – 0795 - May 2023 - Cake - 13-3077	NOT DETECTED	<1.0	-

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm RIDASCREEN FAST Hazelnut R6802	Range for $ z \leq 2$ according to R-Biopharm RIDASCREEN FAST Hazelnut R6802	Result (mg/kg)	Recovery (%)
FAPAS - Hazelnut in Chocolate - 27327A May – July 2022	NOT DETECTED	-	<1.0	-

FAPAS - Hazelnut in Chocolate - 27327B May – July 2022	10.3	5.13 – 15.4	15.3	148.5
FAPAS - Hazelnut and Peanut in Cooked Biscuit -27319 - March – May 2022	18.0	9.01 – 27.0	17.69	98.3
FAPAS - Hazelnut in Chocolate - 27299A - May – July 2021	16.4	8.19 – 24.6	11.22	68.4
FAPAS - Hazelnut in Chocolate - 27299B - May – July 2021	NOT DETECTED	-	<1.0	-
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27291 March – May 2021	10.8	5.42 – 16.3	12.3	113.9

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS - Nuts In Chocolate - 27453 - November 2025 – January 2026	NOT DETECTED	<1.0	97.0
FAPAS - Hazelnut in Chocolate-27433A - May – July 2025	DETECTED	DETECTED	100.0
FAPAS - Hazelnut in Chocolate-27433B - May – July 2025	NOT DETECTED	<1.0	95.0
FAPAS - Hazelnut and Peanut in Cooked Biscuit -27425 - March – April 2025	DETECTED	DETECTED	95.0
BIPEA - Food Allergens - Nuts 77G - 17 - May 2024 - Cake - 12-2877	DETECTED	DETECTED	100.0
BIPEA - Food Allergens - Nuts 77G - 17 - May 2024 - Cake - 10-2977	NOT DETECTED	<1.0	75.0
BIPEA - Food Allergens - Nuts 77G - 17 - May 2024 - Cake - 14-3077	DETECTED	DETECTED	100.0
FAPAS – Hazelnut and Peanut in Cooked Biscuit - 27389 - March – May 2024	DETECTED	5.70	96.0



FAPAS - Nuts In Chocolate - 27378 - November 2023 – January 2024	NOT DETECTED	<1.0	100.0
FAPAS - Nuts In Chocolate - 27343 -November 2022 – January 2023	DETECTED	DETECTED	100.0
FAPAS - Nuts In Chocolate - 27312 -November 2021 – February 2022	NOT DETECTED	<1.0	96.0