

| Proficiency Test | Assigned value (mg/kg) of allergenic protein | Result (mg/kg) | Z-score |
|---|--|----------------|---------|
| FAPAS Lupin in Flour 27388A February – April 2024 | 19.6 | 18.1 | -0.3 |
| FAPAS Lupin in Flour 27388B February – April 2024 | NOT DETECTED | <1.0 | - |

| Proficiency Test | Assigned value (mg/kg) according to R-Biopharm RIDASCREEN FAST Lupine R6102 | Range for $ z \leq 2$ according to R-Biopharm RIDASCREEN FAST Lupine R6102 | Result (mg/kg) | Recovery (%) |
|---|---|---|----------------|--------------|
| FAPAS Lupin in Flour 27424A March – April 2025 | 11.4 | 5.68 – 17.0 | 12.12 | 106.3 |
| FAPAS Lupin in Flour 27424B March – April 2025 | NOT DETECTED | - | <1.0 | - |
| FAPAS Lupin in Flour 27388A February – April 2024 | 20.9 | 10.4 – 31.3 | 18.1 | 86.6 |

| | | | | |
|--|--------------|-------------|-------|-------|
| FAPAS Lupin in Flour 27388B February – April 2024 | NOT DETECTED | - | <1.0 | - |
| FAPAS Lupin in Flour 27352A March – April 2023 | 23.8 | 11.9 – 35.7 | 31.5 | 132.4 |
| FAPAS Lupin in Flour 27352B March – April 2023 | NOT DETECTED | - | <1.0 | - |
| FAPAS Lupin in Flour 27318A March – April 2022 | 19.4 | 9.71 – 29.1 | 18.55 | 95.6 |
| FAPAS Lupin in Flour 27318B March – April 2022 | NOT DETECTED | - | <1.0 | - |