

| Proficiency Test | Assigned value (mg/L) | Result (mg/L) | Z-score |
|--|-----------------------|---------------|---------|
| FAPAS - Hazelnut and Peanut in Cooked Biscuit – 27353 - March – May 2023 | 15.3 | 15.0 | -0.1 |

| Proficiency Test | Assigned value (mg/kg) according to R-Biopharm Ridascreen Peanut R6811 | Range for $ z \leq 2$ according to R-Biopharm Ridascreen Peanut R6811 | Result (mg/kg) | Recovery (%) |
|---|--|--|----------------|--------------|
| FAPAS - Hazelnut and Peanut in Cooked Biscuit -27425 - March – April 2025 | 34.5 | 17.3 – 51.8 | 19.7 | 51.9 |
| FAPAS – Hazelnut and Peanut in Cooked Biscuit - 27389 - March – May 2024 | 7.29 | 3.64 – 10.9 | 10.3 | 141.3 |
| FAPAS Peanut in Garlic Powder 27325A - May – July 2022 | 17.0 | 8.52 – 25.5 | 11.41 | 67.1 |
| FAPAS Peanut in Garlic Powder 27325B - May – July 2022 | NOT DETECTED | - | <1.0 | - |
| FAPAS Hazelnut and Peanut in Cooked Biscuit 27319 - March – May 2022 | 21.0 | 10.5 – 31.5 | 29.71 | 141.5 |
| FAPAS Peanut in Chocolate 27315A - February – March 2022 | NOT DETECTED | - | <1.0 | - |
| FAPAS Peanut in Chocolate 27315B - February – March 2022 | 20.2 | 10.1 – 30.3 | 13.37 | 66.2 |
| FAPAS Peanut in Garlic Powder 27297A - June – August 2021 | 97.0 | 48.5 – 146 | 72.4 | 74.6 |

| | | | | |
|--|--------------|---|------|---|
| FAPAS Peanut in Garlic Powder 27297B - June – August 2021 | NOT DETECTED | - | <1.0 | - |
|--|--------------|---|------|---|

| Proficiency Test | Assignment | Result (mg/kg) | Satisfactory results / results agreeing with assignment (%) |
|---|-------------------|---------------------------|--|
| FAPAS Peanut as the Allergenic Protein in Chocolate 27437A - July – August 2025 | DETECTED | DETECTED | 100.0 |
| FAPAS Peanut as the Allergenic Protein in Chocolate 27437B - July – August 2025 | NOT DETECTED | <0.25 | 100.0 |
| FAPAS Allergens on Swab 27419A - May – July 2025 | DETECTED | DETECTED | 100.0 |
| FAPAS Allergens on Swab 27419B - May – July 2025 | NOT DETECTED | <0.05 | 92.0 |
| FAPAS - Peanut in Chocolate – 27420A - February – March 2025 | NOT DETECTED | <1.0 | 97.0 |
| FAPAS - Peanut in Chocolate – 27420B - February – March 2025 | DETECTED | DETECTED | 97.0 |
| FAPAS Fish and Peanut in Sauce (Cooked) - 27412A – December 2024 – February 2025 | NOT DETECTED | <1.0 | 97.0 |
| FAPAS Fish and Peanut in Sauce (Cooked) – 27412B – December 2024 – February 2025 | DETECTED | DETECTED | 100.0 |
| FAPAS Allergens on Swab 27383A - January – March 2024 | DETECTED | DETECTED | 69.0 |
| FAPAS Allergens on Swab 27383B - January – March 2024 | NOT DETECTED | <1.0 | 94.0 |
| FAPAS Nuts In Chocolate 27312 - November 2021 – February 2022 | DETECTED | DETECTED | 98.0 |