



Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March - April 2026	5.26 (mg/kg of allergenic protein)	5.26	0.0
	18.1	21.0	0.6
FAPAS - Peanut as the Allergenic Protein in Chocolate - 27461A - February - March 2026	NOT DETECTED	<0.25	-
FAPAS - Peanut as the Allergenic Protein in Chocolate - 27461B - February - March 2026	2.93	3.01	0.1
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27353 - March - May 2023	15.3	15.0	-0.1

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Peanut R6811	Range for $z \leq 2$ according to R-Biopharm Ridascreen Peanut R6811	Result (mg/kg)	Recovery (%)
FAPAS - Fish and Peanut in Sauce (Cooked) – 27451A – December 2025 – February 2026	14.9	7.44 – 22.3	14.1	94.6
FAPAS - Fish and Peanut in Sauce (Cooked) – 27451B – December 2025 – February 2026	NOT DETECTED	-	<1.0	-
FAPAS - Hazelnut and Peanut in Cooked Biscuit -27425 - March – April 2025	34.5	17.3 – 51.8	19.7	51.9
FAPAS – Hazelnut and Peanut in Cooked Biscuit - 27389 - March – May 2024	7.29	3.64 – 10.9	10.3	141.3
FAPAS Peanut in Garlic Powder 27325A - May – July 2022	17.0	8.52 – 25.5	11.41	67.1
FAPAS Peanut in Garlic Powder 27325B - May – July 2022	NOT DETECTED	-	<1.0	-
FAPAS Hazelnut and Peanut in Cooked Biscuit 27319 - March – May 2022	21.0	10.5 – 31.5	29.71	141.5
FAPAS Peanut in Chocolate 27315A - February – March 2022	NOT DETECTED	-	<1.0	-
FAPAS Peanut in Chocolate 27315B - February – March 2022	20.2	10.1 – 30.3	13.37	66.2
FAPAS Peanut in Garlic Powder 27297A - June – August 2021	97.0	48.5 – 146	72.4	74.6
FAPAS Peanut in Garlic Powder 27297B - June – August 2021	NOT DETECTED	-	<1.0	-

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Peanut, Sesame & Soya on Swabs (Cotton) 27460A - January - March 2026	NOT DETECTED	<0.1	100.0
FAPAS Peanut, Sesame & Soya on Swabs (Cotton) 27460B - January - March 2026	DETECTED	DETECTED	90.0
FAPAS - Nuts In Chocolate - 27453 - November 2025 - January 2026	NOT DETECTED	<1.0	94.0
FAPAS Peanut as the Allergenic Protein in Chocolate 27437A - July - August 2025	DETECTED	DETECTED	100.0
FAPAS Peanut as the Allergenic Protein in Chocolate 27437B - July - August 2025	NOT DETECTED	<1.0	100.0
FAPAS Allergens on Swab 27419A - May - July 2025	DETECTED	DETECTED	100.0
FAPAS Allergens on Swab 27419B - May - July 2025	NOT DETECTED	<0.05	92.0
FAPAS - Peanut in Chocolate - 27420A - February - March 2025	NOT DETECTED	<1.0	97.0
FAPAS - Peanut in Chocolate - 27420B - February - March 2025	DETECTED	DETECTED	97.0
FAPAS Fish and Peanut in Sauce (Cooked) - 27412A - December 2024 - February 2025	NOT DETECTED	<1.0	97.0
FAPAS Fish and Peanut in Sauce (Cooked) - 27412B - December 2024 - February 2025	DETECTED	DETECTED	100.0
FAPAS Allergens on Swab 27383A - January - March 2024	DETECTED	DETECTED	69.0
FAPAS Allergens on Swab 27383B - January - March 2024	NOT DETECTED	<1.0	94.0
FAPAS Nuts In Chocolate 27312 - November 2021 - February 2022	DETECTED	DETECTED	98.0

