


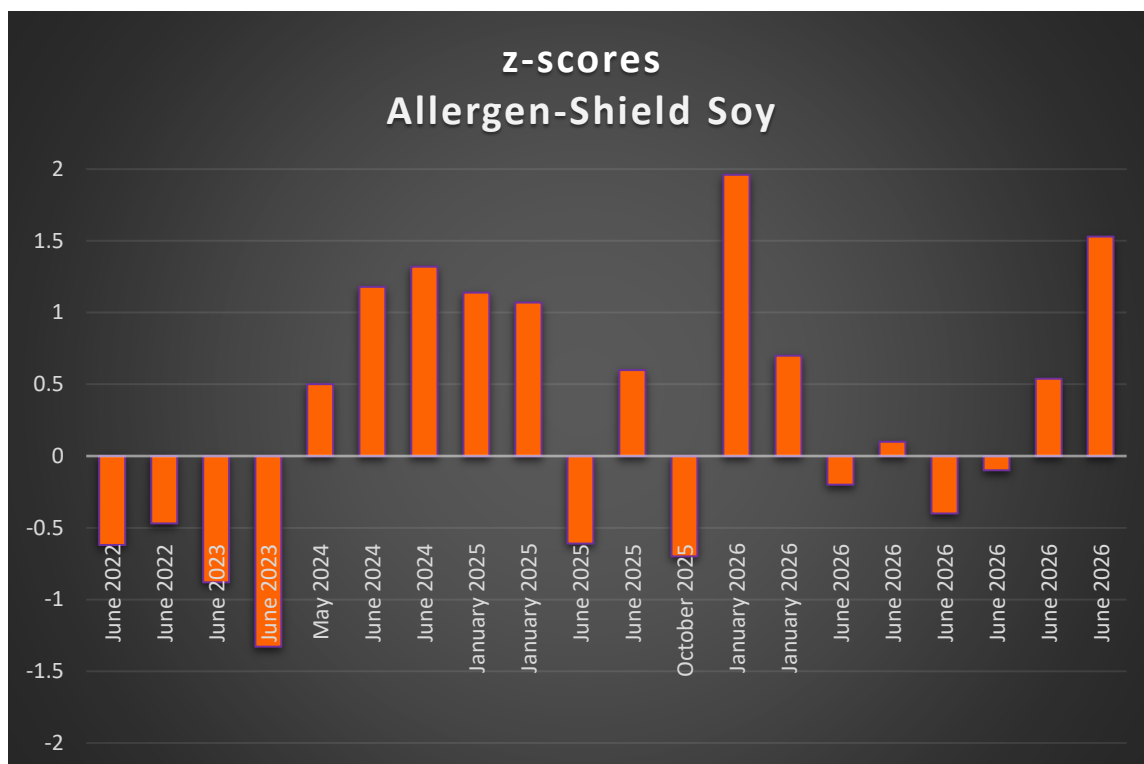

Proficiency Test	Assigned value (mg/kg)	Result (mg/kg)	Z-score
BIPEA Food Allergens - Soya - 77H - Wheat Flour - 10 - 3177 - June 2026	31.0	35.9	0.54
BIPEA Food Allergens - Soya - 77H - Wheat Flour - 10 - 3277 - June 2026	16.4	22.9	1.53
FAPAS Soya in Wheat Flour 27469A - April - May 2026	30.1	28.5	-0.2
FAPAS Soya in Wheat Flour 27469A - April - May 2026	27.9	28.5	0.1
FAPAS Soya in Wheat Flour 27469B - April - May 2026	22.7 (allergenic protein)	20.5	-0.4

FAPAS Soya in Wheat Flour 27469B - April – May 2026	21.2	20.5	-0.1
BIPEA Food Allergens - Soya - 77H – Meat - 11 - 2277 - January 2026	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Soya - 77H – Meat - 11 - 2377 - January 2026	6.40	10.8	1.96
BIPEA Food Allergens - Soya - 77H – Meat - 11 - 2477 - January 2026	32.7	39.8	0.70
FAPAS Soya in Wheat Flour 27446A - September – October 2025	NOT DETECTED	<1.0	-
FAPAS Soya in Wheat Flour 27446B - September – October 2025	30.3	24.7	-0.7
BIPEA Food Allergens - Soya - 77H – Wheat Flour - 09 - 3177 - June 2025	50.4	44.6	-0.61
BIPEA Food Allergens - Soya - 77H – Wheat Flour - 09 - 3277 - June 2025	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Soya - 77H – Wheat Flour - 10 - 3377 - June 2025	18.3	19.2	0.60
BIPEA Food Allergens - Soya - 77H – Meat - 10 - 2277 - January 2025	58.4	82.4	1.14
BIPEA Food Allergens - Soya - 77H – Meat - 10 - 2377 - January 2025	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Soya - 77H – Meat - 10 - 2477 - January 2025	27.2	38.4	1.07
BIPEA Food Allergens - Soya - 77H – Wheat Flour - 09 - 3377 - June 2024	39.1	54.1	1.32
BIPEA Food Allergens - Soya - 77H – Wheat Flour - 08 - 3277 - June 2024	NOT DETECTED	<2.5	-

BIPEA Food Allergens - Soya - 77H - Wheat Flour - 08 - 3177 - June 2024	14.1	18.7	1.18
FAPAS Soya in Wheat Flour 27392B - April - May 2024	21.9	24.5	0.5
BIPEA Food Allergens - Soya - 77H - Flour - 07 - 3177 - June 2023	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Soya - 77H - Flour - 07 - 3277 - June 2023	64.0	42.8	-1.33
BIPEA Food Allergens - Soya - 77H - Flour - 08 - 3377 - June 2023	27.5	20.7	-0.88
BIPEA Food Allergens - Soya - 77H - Flour - 06 - 3177 - June 2022	15.7	13.2	-0.47
BIPEA Food Allergens - Soya - 77H - Flour - 06 - 3277 - June 2022	37.3	32.1	-0.62
BIPEA Food Allergens - Soya - 77H - Flour - 06 - 3377 - June 2022	NOT DETECTED	<2.5	-

Proficiency Test	Assigned value (mg/kg) according to R- Biopharm Ridascreen Fast Soya R7102	Range for $ z \leq 2$ according to R- Biopharm Ridascreen Fast Soya R7102	Result (mg/kg)	Recovery (%)
FAPAS Soya in Wheat Flour 27356A - May - June 2023	NOT DETECTED (89% CONSENSUS)	-	<1.0	-
FAPAS Soya in Wheat Flour 27356B - May - June 2023	29.0	14.5 - 43.5	22.5	77.6
FAPAS Soya in Wheat Flour 27337A - September - November 2022	DETECTED (VALUE NOT SET)	-	9.18	-

FAPAS Soya in Wheat Flour 27337B - September - November 2022	20.9	10.4 – 31.3	17.38	83.2
FAPAS Soya in Wheat Flour 27323A - April – June 2022	DETECTED	-	DETECTED	-
FAPAS Soya in Wheat Flour 27323B - April – June 2022	15.2	7.58 – 22.8	21.0	138.2
FAPAS Soya in Wheat Flour 27295A - April – May 2021	10.4	5.19 – 15.6	11.49	110.5
FAPAS Soya in Wheat Flour 27295B - April – May 2021	26.0	13.0 – 39.0	26.32	101.2

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Peanut, Sesame & Soya on Swabs (Cotton) 27460E - January – March 2026	NOT DETECTED	<0.1	90.0
FAPAS Peanut, Sesame & Soya on Swabs (Cotton) 27460F - January – March 2026	DETECTED	DETECTED	45.0
FAPAS Allergens on Swab 27419E - May – July 2025	NOT DETECTED	<0.05	91.0
FAPAS Allergens on Swab 27419F - May – July 2025	DETECTED	DETECTED	39.0
FAPAS Soya in Wheat Flour 27392A - April – May 2024	DETECTED	DETECTED	100.0
FAPAS Soya in Wheat Flour 27392B - April – May 2024	DETECTED	DETECTED	100.0

FAPAS Allergens on Swab 27383E - January – March 2024	NOT DETECTED	NOT DETECTED	100.0
FAPAS Allergens on Swab 27383F - January – March 2024	DETECTED	DETECTED	63.0
BIPEA Food Allergens - Soya - 77H - Meat - 09 - 2277 – January 2024	DETECTED	DETECTED	100.0
BIPEA Food Allergens - Soya - 77H - Meat - 09 - 2377 – January 2024	DETECTED	DETECTED	100.0
BIPEA Food Allergens - Soya - 77H - Meat - 10 - 2477 – January 2024	NOT DETECTED	<0.2	100.0
BIPEA Food Allergens - Soya - 77H - Meat - 08 - 2277 – January 2023	DETECTED	DETECTED	100.0
BIPEA Food Allergens - Soya - 77H - Meat - 08 - 2377 – January 2023	NOT DETECTED	<0.2	100.0
BIPEA Food Allergens - Soya - 77H - Meat - 09 - 2477 – January 2023	DETECTED	DETECTED	100.0