



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31E - Mycotoxins - Feed - Animal Feed - 14 - 5 - January 2024	5.83	4.68	-0.67
BIPEA 31B - Mycotoxins - Cereals - Ground Corn - 6 - 187 - December 2023	5.87	8.78	1.69
FAPAS Food Chemistry Proficiency Test 04497 Aflatoxins in Peanut Butter (Water Slurry) November 2023 - January 2024	3.58	3.89	0.4



BIPEA 31A - Mycotoxins - Dried Fruits, Spices and Other Products - Peanut Paste - 6 - 185 - September 2023	8.69	9.57	0.36
FAPAS Food Chemistry Proficiency Test 04482 Multi-Mycotoxins in Maize Flour July - September 2023	4.72	4.73	0.0
FAPAS Food Chemistry Proficiency Test 04472 Multi-Mycotoxins in Maize Flour April - June 2023	4.73	4.39	-0.3
FAPAS Food Chemistry Proficiency Test 22199 Multi-Mycotoxins in Wheat Flour November 2022-January 2023	3.78	3.34	-0.5
FAPAS Food Chemistry Proficiency Test 04449 Mycotoxins in Dried Figs (Water/Fruit Slurry) August - September 2022	3.32	3.52	0.3
FAPAS Food Chemistry Proficiency Test 04448 Multi-Mycotoxins in Maize Flour July - September 2022	3.32	3.71	0.5
FAPAS Food Chemistry Proficiency Test 04439 Multi-Mycotoxins in Maize Flour April - May 2022	5.55	5.60	0.0
FAPAS Food Chemistry Proficiency Test 04438 Aflatoxins in Animal Feed (Cereal Based) April - May 2022	7.80	7.75	0.0



	Negative Sample	<limit of detection	-
A.I.A. Ring Test Lotto RTB1 220322 Aflatoxin B1 in Corn March 2022	9.22	10.17	0.32
	17.76	19.26	0.31
			0.52
FAPAS Food Chemistry Proficiency Test 22187 Multi-Mycotoxins in Wheat Flour November 2021 - January 2022	2.60	2.88	0.5
FAPAS Food Chemistry Proficiency Test 04426 Multi-Mycotoxins in Maize Flour October - December 2021	5.82	6.37	0.3
FAPAS Food Chemistry Proficiency Test 04418 Multi-Mycotoxins in Maize Flour July - September 2021	3.12	3.35	0.3
FAPAS Food Chemistry Proficiency Test 04417 Mycotoxins in Hazelnut (Water/Nut Slurry) June - August 2021	4.20	4.81	0.7
FAPAS Food Chemistry Proficiency Test 04411 Multi-Mycotoxins in Maize Flour April - June 2021	5.35	5.25	-0.1
FAPAS Food Chemistry Proficiency Test 04408 Aflatoxins in Peanut	4.86	5.64	0.7



(Water/Nut Slurry) March - April 2021			
FAPAS Food Chemistry Proficiency Test 04402 Mycotoxin Contamination in Animal Feed (Cereal Based) December 2020 - January 2021	20.4	18.18	-0.5
FAPAS Food Chemistry Proficiency Test 04390 Mycotoxins in Hazelnut (Water/Nut Slurry) July - August 2020	4.50	3.67	-0.8
FAPAS Food Chemistry Proficiency Test 04384 Multi-Mycotoxins in Maize Flour April - May 2020	6.13	5.95	-0.1
FAPAS Food Chemistry Proficiency Test 04378 Multi-Mycotoxins in Maize Flour February - March 2020	6.65	6.48	-0.1
FAPAS Food Chemistry Proficiency Test 04375 Mycotoxin Contamination in Animal Feed (Cereal Based) December 2019 - January 2020	14.6	12.78	-0.6
A.I.A. Ring Test Lotto RTB1 240919 Aflatoxin B1 in Corn September 2019	<limit of detection	<limit of detection	- 0.142
	7.2	7.15	
	13.9	12.90	
	22.3	22.07	

