



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31A – 5 - 187 - Mycotoxins – Dried Fruits, Spices and Other Products – Coffee – 5 - 1731 – January 2024	7.89	6.38	-0.68
FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 – January 2024	2.31	2.23	-0.2
FAPAS Food Chemistry Proficiency Test - 22215 - Multi - Mycotoxins in Wheat Flour - November 2023 – January 2024	1.91	2.27	0.9



FAPAS Food Chemistry Proficiency Test - 04496 - Aflatoxins in Rice – November 2023 – January 2024	4.13	3.19	-1.0
BIPEA 31E - 186 - Mycotoxins - Feed – Corn Silage – 15 - 5831 - November 2023	5.48	4.64	-0.52
BIPEA 31A - 186 - Mycotoxins - Dried Fruits, Spices and Other Products – Ground Ginger – 02 - 1231 - November 2023	7.97	8.00	0.01
BIPEA 31B – 186 - Mycotoxin Contamination in Cereals – Common Wheat – 3 - 0631 - October 2023	3.30	2.63	-0.68
FAPAS Food Chemistry Proficiency Test - 04481 - Mycotoxins in Pet Dog Food - July – August 2023	9.88	5.66	-1.9
BIPEA 31B – 185 - Mycotoxin Contamination in Cereals – Rice - 06-4331 - June 2023	4.47	3.14	-0.99
BIPEA 31A - 184 Mycotoxins - Dried Fruits, Spices and Other Products – Pistachio Paste – 08 - 1431 - May 2023	10.05	13.45	1.23
A.I.A. Ring Test Lotto - RT B1 300523 - Aflatoxin B1 in Corn - May 2023	3.51	2.76	-0.57
	7.92	12.46	1.57
	15.08	5.86	-1.80
	27.21	26.00	-0.16



BIO-SHIELD

BIPEA 31B - 184 - Mycotoxin Contamination in Cereals - Rusk (Crispbread) 02-5431 - April 2023		2.65	3.63	1.23
BIPEA 31E - 3 - Mycotoxin Contamination in Animal feed - 13 - 3031 - January 2023		3.16	3.65	0.52
BIPEA 31B - 183 - Mycotoxin Contamination in Ground Barley - 05 - 0531 - December 2022		6.42	6.62	0.11
BIPEA 31B - 182 - Mycotoxin Contamination in Ground Corn - 18 - 0731 - December 2022		5.7	2.9	-1.65
BIPEA 31B - 180 - Mycotoxin Contamination in Rice - 05 - 4331 - June 2022		2.64	3.65	1.28
BIPEA 31E - 2 - Mycotoxin Contamination in Corn Silage - 01 - 5831 - May 2022		2.16	1.12	-1.6
FAPAS Food Chemistry Proficiency Test - 04430 - Mycotoxin Contamination in Animal Feed - December 2021 - January 2022		19.1	19.5	0.5
QUALINOVA ROUND : AL 21008 - Aflatoxins in dried fruit - April 2021		3.95	3.29	-0.76
DRRR - RVEP 190663 - CORN AFLATOXIN 1	Sample 1	8.6	9.4	0.53
	Sample 2	15.2	15.1	-0.03



FAPAS Food Chemistry Proficiency Test - 04366 - Multi - Mycotoxins in Maize - August - September 2019	1.88	1.84	-0.1
FAPAS Food Chemistry Proficiency Test - 04359 - Multi - Mycotoxins in Maize - April - May 2019	9.22	9.58	0.2
A.I.A. Ring Test Lotto - RT B1 260319 - Aflatoxin B1 in Corn - April 2019	<limit of detection	<0.5	0.117
	7.60	8.94	
	14.10	15.69	
	29.40	28.11	
FAPAS Food Chemistry Proficiency Test - 04354 - Multi - Mycotoxins in Maize - February - March 2019	4.85	4.58	-0.3
FAPAS Food Chemistry Proficiency Test - 04330 - Multi - Mycotoxins in Maize - February - March 2018	3.83	3.39	-0.5
FAPAS Food Chemistry Proficiency Test - 04319 - Multi - Mycotoxins in Maize - August - September 2017	4.69	5.12	0.4