



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 - January 2024	1240	1224	-0.1
FAPAS Food Chemistry Proficiency Test - 04498 - Mycotoxin Contamination in Animal Feed (Cereal Based) - November 2023 - January 2024	2800	2759	-0.1
FAPAS Food Chemistry Proficiency Test - 22215 - Multi - Mycotoxins in Wheat Flour - November 2023 - January 2024	830	872.9	0.3



BIO-SHIELD

BIPEA 31B - Mycotoxins - Cereals - Ground Corn - 6 - 187 - December 2023	1918	1464	-1.11
FAPAS Food Chemistry Proficiency Test - 04482 - Multi - Mycotoxins in Maize Flour July - September 2023	1380	1238	-0.7
BIPEA 31E - 3 Mycotoxin Contamination in Animal feed - 13-3031 - January 2023	939	975	0.17
FAPAS Food Chemistry Proficiency Test - 04461 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2022 - January 2023	Not set	<100	-
FAPAS Food Chemistry Proficiency Test - 22199 - Multi - Mycotoxins in Wheat Flour - November 2022-January 2023	809	732.50	-0.6
FAPAS Food Chemistry Proficiency Test - 22187 - Multi - Mycotoxins in Wheat Flour November 2021 - January 2022	745	768.58	0.2
FAPAS Food Chemistry Proficiency Test - 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021	959	1088	0.8
FAPAS Food Chemistry Proficiency Test - 22177 - Multi - mycotoxins in Wheat Flour - November 2020 - January 2021	769	731.20	-0.3
FAPAS Food Chemistry Proficiency Test - 22175 - Deoxynivalenol, Zearalenone, T-2 and HT-2 in Animal Feed - September - October 2020	900	877.52	-0.2



BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04384 - Multi - Mycotoxins in Maize Flour - April - May 2020	859	847.8	-0.1
FAPAS Food Chemistry Proficiency Test - 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	1200	1218.22	0.1
FAPAS Food Chemistry Proficiency Test - 04375 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2019 - January 2020	2702	2320.55	-1.0
FAPAS Food Chemistry Proficiency Test - 22166 - Deoxynivalenol, Zearalenone, T-2 and HT-2 in Wheat Flour - November 2019 - January 2020	708	654.28	-0.4