


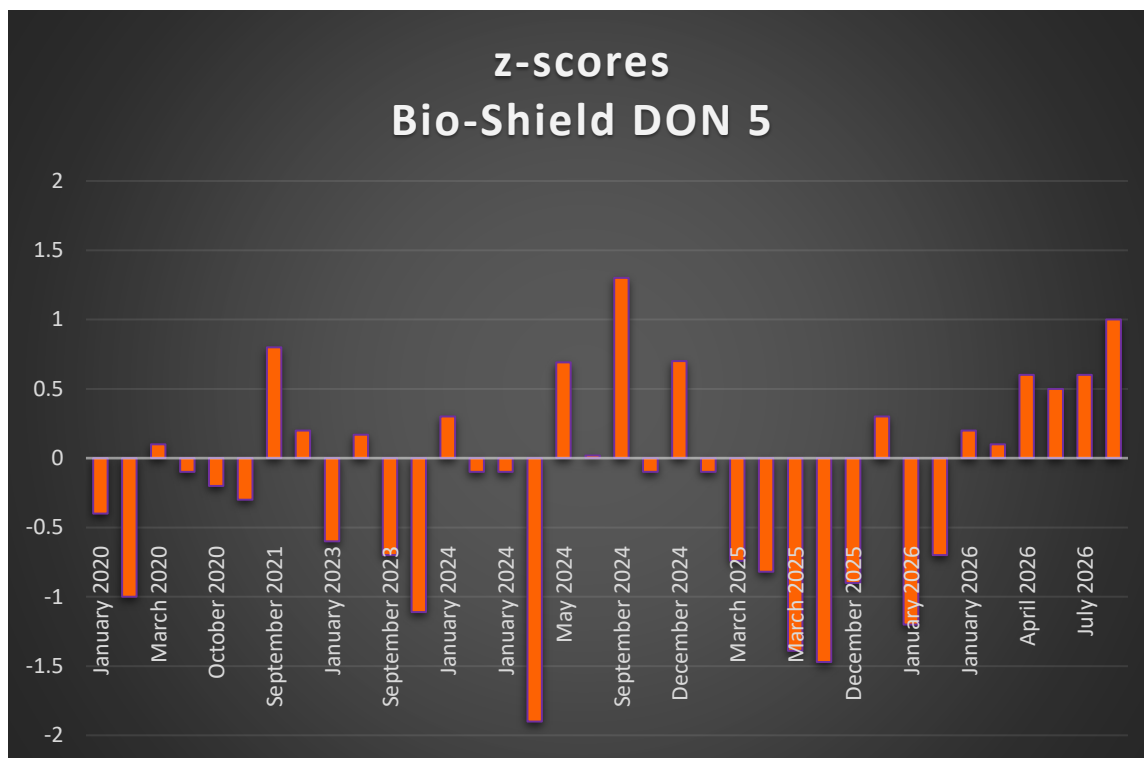

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2026-R01-Tx01] [Wheat Flour] – July 2026	11.6	34.75	0.5
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2026-R01-Tx02] [Wheat Flour] – July 2026	490.7	569.95	0.6
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2026-R01-Tx03] [Wheat Flour] – July 2026	794.6	995.51	1.0
BIPEA 31B – Mycotoxins - Cereals – Pasta - 1 - 6431 – April 2026	202	233	0.6



FAPAS Food Chemistry Proficiency Test - 04578 - Multi-mycotoxins in Durum Wheat - November 2025 - January 2026	825	841	0.1
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R07-Tx01] [Wheat] - January 2026	1132.75	903	-1.2
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R07-Tx02] [Wheat] - January 2026	519	440	-0.7
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R07-Tx03] [Wheat] - January 2026	18.2	27.7	0.2
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R04-Tx01] [Wheat] - December 2025	827	696	-0.9
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R04-Tx02] [Wheat] - December 2025	195	211	0.3
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R04-Tx03] [Wheat] - December 2025	13.0	<100	-
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx01] [Barley] - March 2025	396.8	323	-0.74
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx02] [Wheat] - March 2025	1223.4	1059	-0.82



UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx03] [Wheat] – March 2025	11.0	<100	-
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx01] [Wheat] – March 2025	1046	795.2	-1.39
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx02] [Wheat] – March 2025	1046	780.9	-1.47
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx03] [Wheat] – March 2025	15.0	<100	-
FAPAS Food Chemistry Proficiency Test - 04540 – Mycotoxins in Maize Flour - January – March 2025	750	737	-0.1
BIPEA 31C - Mycotoxins – Baby Food –24- 3931 – December 2024	159	188	0.70
FAPAS Food Chemistry Proficiency Test - 04519 – Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July - September 2024	803	790	-0.1
FAPAS Food Chemistry Proficiency Test - 04520 - Multi - Mycotoxins in Maize Flour - July – September 2024	782	951	1.3
BIPEA 31C - Mycotoxins – Baby Food –23 - 3931 – June 2024	204	205	0.02
BIPEA 31E - Mycotoxins – Feed – Soyabean Meal - 16 - 6031 – May 2024	219	258	0.69



FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April – May 2024	759	523.4	-1.9
FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 – January 2024	1240	1224	-0.1
FAPAS Food Chemistry Proficiency Test - 04498 – Mycotoxin Contamination in Animal Feed (Cereal Based) - November 2023 – January 2024	2800	2759	-0.1
FAPAS Food Chemistry Proficiency Test - 22215 - Multi - Mycotoxins in Wheat Flour - November 2023 – January 2024	830	872.9	0.3
BIPEA 31B - Mycotoxins – Cereals – Ground Corn - 6 - 187 – December 2023	1918	1464	-1.11
FAPAS Food Chemistry Proficiency Test - 04482 - Multi - Mycotoxins in Maize Flour July - September 2023	1380	1238	-0.7
BIPEA 31E - 3 Mycotoxin Contamination in Animal feed - 13-3031 - January 2023	939	975	0.17
FAPAS Food Chemistry Proficiency Test - 04461 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2022 - January 2023	Not set	<100	-
FAPAS Food Chemistry Proficiency Test - 22199 – Multi - Mycotoxins in Wheat Flour - November 2022-January 2023	809	732.50	-0.6



FAPAS Food Chemistry Proficiency Test - 22187 - Multi - Mycotoxins in Wheat Flour November 2021 – January 2022	745	768.58	0.2
FAPAS Food Chemistry Proficiency Test - 04418 - Multi – Mycotoxins in Maize Flour - July – September 2021	959	1088	0.8
FAPAS Food Chemistry Proficiency Test - 22177 - Multi - mycotoxins in Wheat Flour - November 2020 - January 2021	769	731.20	-0.3
FAPAS Food Chemistry Proficiency Test - 22175 - Deoxynivalenol, Zearalenone, T-2 and HT-2 in Animal Feed - September – October 2020	900	877.52	-0.2
FAPAS Food Chemistry Proficiency Test – 04384 - Multi – Mycotoxins in Maize Flour - April – May 2020	859	847.8	-0.1
FAPAS Food Chemistry Proficiency Test – 04378 - Multi – Mycotoxins in Maize Flour - February – March 2020	1200	1218.22	0.1
FAPAS Food Chemistry Proficiency Test - 04375 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2019 - January 2020	2702	2320.55	-1.0
FAPAS Food Chemistry Proficiency Test - 22166 - Deoxynivalenol, Zearalenone, T-2 and HT-2 in Wheat Flour - November 2019 - January 2020	708	654.28	-0.4