



Test	Assigned value (mg/kg)	Result (mg/kg)	Z-score
BIPEA Food Allergens - Histamine - 77E - Food Allergens - Histamine in Fish 17-1377 - March 2024	94.5	85.7	-0.63
BIPEA Food Allergens - Histamine - 77E - Food Allergens - Histamine in Fish 18-1477 - March 2024	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens - Histamine in Fish 19-1577 - March 2024	28.7	30.4	0.16
Lab Conformance Dosage De l'histamine et de l'ABVT dans les produits de la	33.0	34.5	0.4



BIO-SHIELD

peche – AR-23 B - Niveau 1 - Octobre 2023			
Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – AR-23 B - Niveau 2 - Octobre 2023	81.0	85.7	1.1
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 16-1377 - October 2023	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 17-1477 - October 2023	61.5	57.8	-0.27
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 18-1577 - October 2023	21.4	22.0	0.14
FAPAS Food Chemistry Proficiency Test 27367 Histamine in Canned Fish - August – September 2023	27.6	26.2	-0.5
Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – Niveau 1 - Juin 2023	22.7	21.4	-0.3
Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – Niveau 2 - Juin 2023	59.3	61.0	0.2
FAPAS Food Chemistry Proficiency Test 27354 Histamine in Canned Fish - March – May 2023	120	125.1	0.6



BIO-SHIELD

BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 15- 1377 - March 2023	NOT DETECTED	<2.0	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 16- 1477 - March 2023	45.6	18.9	-1.91
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 17- 1577 - March 2023	19.9	10.3	-1.64
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 14- 1377 - October 2022	24.4	23.7	-0.28
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 15- 1477 - October 2022	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 16- 1577 - October 2022	69.2	70.0	0.06
FAPAS Food Chemistry Proficiency Test 27320 Histamine in Canned Fish - March – May 2022	168	162.6	-0.4
FAPAS Food Chemistry Proficiency Test 27303 Histamine in Canned Fish - August – September 2021	31.3	32.04	0.2
FAPAS Food Chemistry Proficiency Test 27292 Histamine in Canned Fish - March – May 2021	268	258.9	-0.5



BIO-SHIELD

FAPAS Food Chemistry Proficiency Test 27267 Histamine in Canned Fish - March - May 2020	294	317.7	1.2
FAPAS Food Chemistry Proficiency Test 27253 Histamine in Canned Fish - August - September 2019	38.5	39.53	0.30