


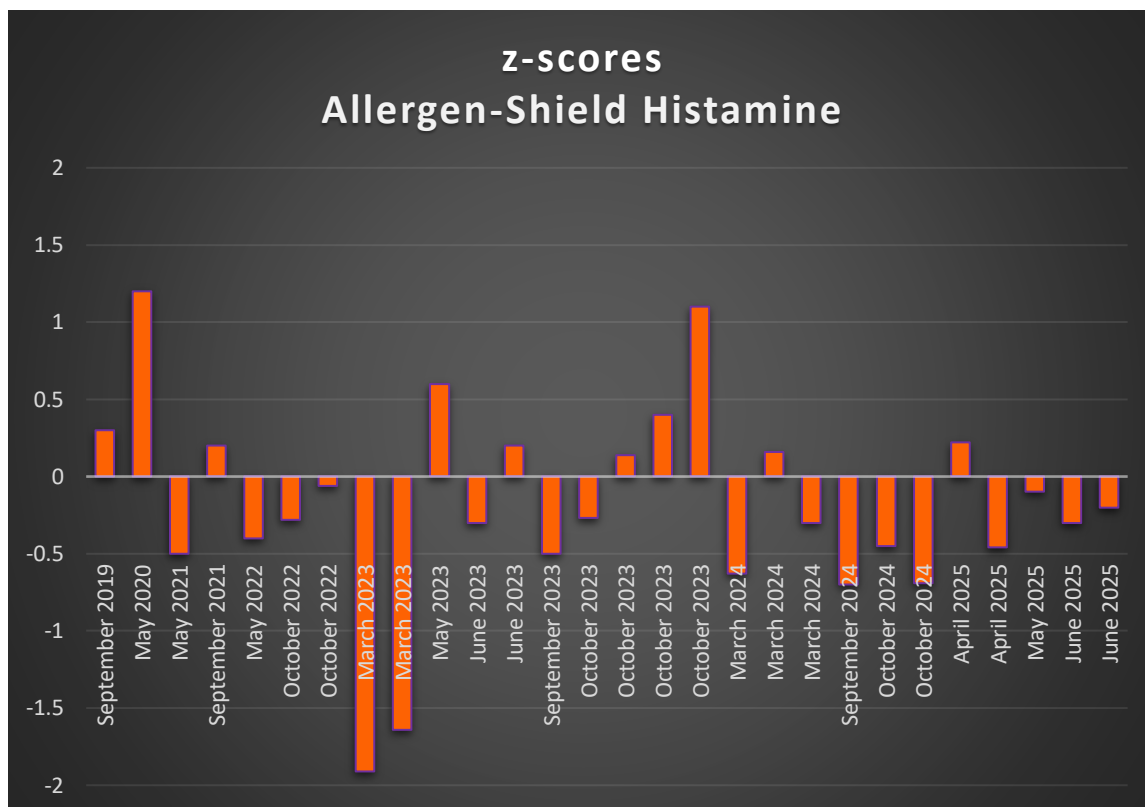

Test	Assigned value (mg/kg)	Result (mg/kg)	Z-score
Lab Conformance Dosage De l'histamine et de l'ABVT dans les produits de la pêche - AR-25 A - Niveau 1 - Juin 2025	22.5	21.4	-0.3
Lab Conformance Dosage De l'histamine et de l'ABVT dans les produits de la pêche - AR-25 A - Niveau 2 - Juin 2025	75.1	74.5	-0.2
FAPAS Food Chemistry Proficiency Test 27426 Histamine in Canned Fish - March - May 2025	81.8	81.2	-0.1



BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 21-1577 – March 2025	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 20-1477 – March 2025	22.3	20.0	-0.46
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 19-1377 – March 2025	51.4	53.8	0.22
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 20-1577 - October 2024	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 19-1477 - October 2024	45.2	35.5	-0.69
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 18-1377 - October 2024	25.8	21.6	-0.45
FAPAS Food Chemistry Proficiency Test 27403 Histamine in Canned Fish - August – September 2024	36.0	33.5	-0.7
FAPAS Food Chemistry Proficiency Test 27390 Histamine in Canned Fish - March – May 2024	273	267	-0.3
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 17-1377 - March 2024	94.5	85.7	-0.63



BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 18-1477 - March 2024	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 19-1577 - March 2024	28.7	30.4	0.16
Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – AR-23 B - Niveau 1 - Octobre 2023	33.0	34.5	0.4
Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – AR-23 B - Niveau 2 - Octobre 2023	81.0	85.7	1.1
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 16-1377 - October 2023	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 17-1477 - October 2023	61.5	57.8	-0.27
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 18-1577 - October 2023	21.4	22.0	0.14
FAPAS Food Chemistry Proficiency Test 27367 Histamine in Canned Fish - August – September 2023	27.6	26.2	-0.5
Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – Niveau 1 - Juin 2023	22.7	21.4	-0.3



Lab Conformite Dosage De l'histamine et de l'ABVT dans les produits de la peche – Niveau 2 - Juin 2023	59.3	61.0	0.2
FAPAS Food Chemistry Proficiency Test 27354 Histamine in Canned Fish - March – May 2023	120	125.1	0.6
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 15-1377 - March 2023	NOT DETECTED	<2.0	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 16-1477 - March 2023	45.6	18.9	-1.91
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 17-1577 - March 2023	19.9	10.3	-1.64
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 14-1377 - October 2022	24.4	23.7	-0.28
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 15-1477 - October 2022	NOT DETECTED	<2.5	-
BIPEA Food Allergens - Histamine - 77E - Food Allergens – Histamine in Fish 16-1577 - October 2022	69.2	70.0	0.06
FAPAS Food Chemistry Proficiency Test 27320 Histamine in Canned Fish - March – May 2022	168	162.6	-0.4



FAPAS Food Chemistry Proficiency Test 27303 Histamine in Canned Fish - August – September 2021	31.3	32.04	0.2
FAPAS Food Chemistry Proficiency Test 27292 Histamine in Canned Fish - March – May 2021	268	258.9	-0.5
FAPAS Food Chemistry Proficiency Test 27267 Histamine in Canned Fish - March – May 2020	294	317.7	1.2
FAPAS Food Chemistry Proficiency Test 27253 Histamine in Canned Fish - August – September 2019	38.5	39.53	0.30