



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31A - 188 - Mycotoxins - Dried Fruits, Spices and Other Products - Hazelnut Paste - 3 - 2031 - March 2024	11.20	8.81	-0.78
BIPEA 31A - 5 - 187 - Mycotoxins - Dried Fruits, Spices and Other Products - Coffee - 5 - 1731 - January 2024	13.24	9.60	-1.02
FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 - January 2024	2.72	3.52	1.3
BIPEA 31C - Mycotoxins - Baby Food - 22 - 22 - 3931 - December 2023	1.62	1.60	-0.04



BIO-SHIELD

BIPEA 31E - 186 - Mycotoxins - Feed – Corn Silage – 15 - 5831 - November 2023	5.73	7.29	0.93
BIPEA 31A - 186 - Mycotoxins - Dried Fruits, Spices and Other Products – Ground Ginger – 02 - 1231 - November 2023	8.04	9.36	0.58
BIPEA 31B - 186 - Mycotoxin Contamination in Cereals – Common Wheat – 3 - 0631 - October 2023	5.25	5.61	0.23
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Peanut Paste - 6 - 185 – September 2023	10.86	9.57	-0.44
FAPAS Food Chemistry Proficiency Test - 04472 – Multi - Mycotoxins in Maize - April - June 2023	4.12	3.84	-0.3
BIPEA 31B - 184 - Mycotoxin Contamination in Cereals 02 - 5431 - April 2023	3.73	4.38	0.58
BIPEA 31A - 182 - Mycotoxin Contamination in Raisins 07 - 3131 - January 2023	14.7	8.7	-1.52
BIPEA 31C - 20 - Mycotoxins - Baby Food - 20 - 3931 – December 2022	2.30	1.44	-1.25
BIPEA 31B - 182 - Mycotoxin Contamination in Ground corn – 18 - 0731 - December 2022	5.98	3.22	-1.58
BIPEA 31B - 180 - Mycotoxin Contamination in Rice – 05 - 4331 - June 2022	6.76	4.78	-1.02



BIPEA 31E - 2 - Mycotoxin Contamination in Corn Silage - 01-5831 - May 2022	6.72	4.83	-0.98
FAPAS Food Chemistry Proficiency Test - 04430 - Mycotoxin Contamination in Animal Feed - December 2021 - January 2022	74.50	66.85	-0.5
FAPAS Food Chemistry Proficiency Test - 04367 - Mycotoxins in Dried Figs - August - October 2019	4.75	5.23	0.5
FAPAS Food Chemistry Proficiency Test - 04366 - Multi - Mycotoxins in Maize - August - September 2019	1.99	2.04	0.1
FAPAS Food Chemistry Proficiency Test - 04364 - Mycotoxins in Paprika - June - August 2019	22.3	20.2	-0.4
FAPAS Food Chemistry Proficiency Test - 04359 - Multi - Mycotoxins in Maize - April - May 2019	2.69	2.74	0.1
FAPAS Food Chemistry Proficiency Test - 17192 - Ochratoxin A in Dried Vine Fruit (Water/Fruit Slurry) - April - May 2019	8.38	8.27	-0.1
FAPAS Food Chemistry Proficiency Test - 04354 - Multi - Mycotoxins in Maize - February - March 2019	4.43	4.88	0.5
FAPAS Food Chemistry Proficiency Test - 04351 - Mycotoxin Contamination in Animal Feed (cereal based) - December 2018 - January 2019	30.2	28.2	-0.3



BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04330 - Multi - Mycotoxins in Maize - February - March 2018	2.75	2.75	0.0
FAPAS Food Chemistry Proficiency Test - 04319 - Multi-Mycotoxins in Maize - August - September 2017	4.73	5.40	0.6
BIPEA 31B - 154 - Mycotoxins - Cereals - Bran - 02-3631 - April 2017	9.60	8.90	-0.26
BIPEA 31B - 153 - Mycotoxins - Cereals - Animal Feed - 09-3031 - February 2017	8.70	9.03	0.24
FAPAS Food Chemistry Proficiency Test - 04301 - Multi Mycotoxins in Maize - November - December 2016	3.03	3.28	0.4