


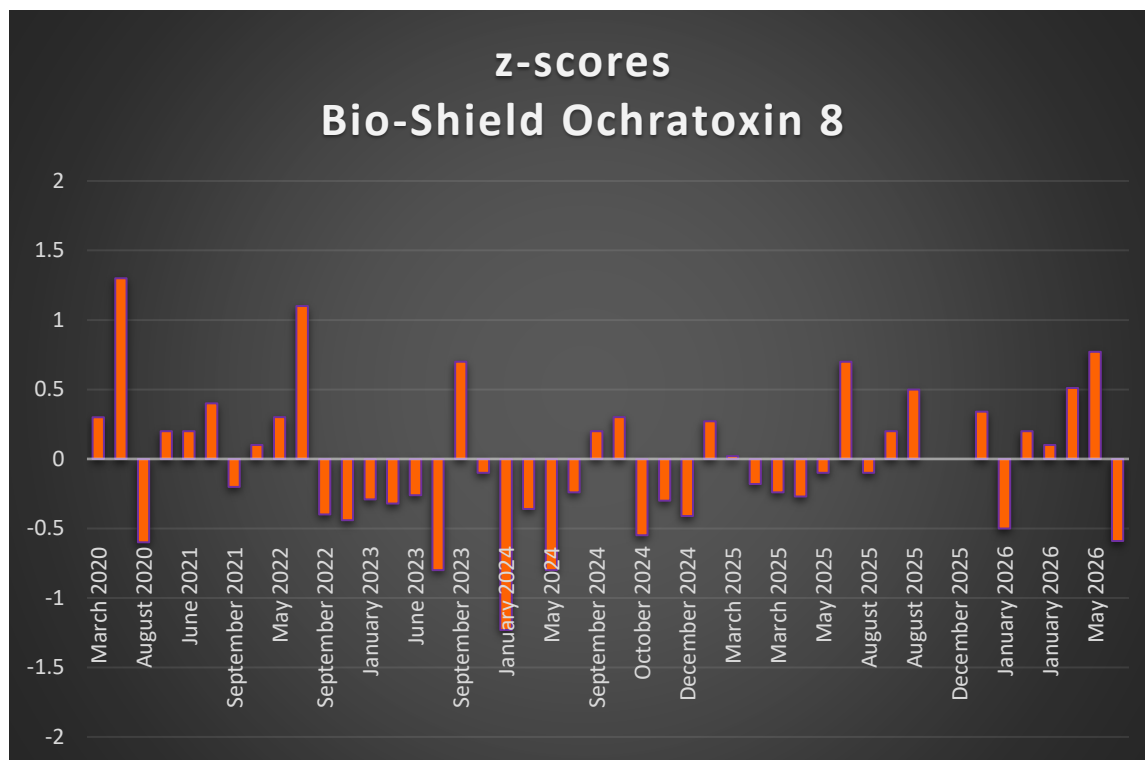

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31B - Mycotoxins – Cereals – Rice – 7 - 4331 – June 2026	7.05	5.87	-0.59
BIPEA 31E - Mycotoxins – Feed – Soyabean Meal – 18 - 6031 – May 2026	5.07	5.84	0.51
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Lentils – 03 - 5931 – May 2026	4.35	5.35	0.77
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R07-Tx01] [Wheat] – January 2026	9.62	8.2	-0.5



UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R07-Tx02] [Wheat] – January 2026	7.013	7.58	0.2
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R07-Tx03] [Wheat] – January 2026	4.555	4.68	0.1
BIPEA 31B - Mycotoxins – Cereals – Ground Corn – 19 - 0731 – December 2025	5.58	6.14	0.34
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R04-Tx01] [Wheat] – December 2025	7.89	8.0	0.0
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R04-Tx02] [Wheat] – December 2025	13.6	13.4	0.0
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2025-R04-Tx03] [Wheat] – December 2025	0.08	<2.0	-
FAPAS Food Chemistry Proficiency Test - 04559 – Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July – August 2025	9.33	10.4	0.5
UK NIR Grain Network Monthly Check Sample Results - HY2025-R01 – Wheat Flour (Spiked) – August 2025	0.69	1.05	0.7
UK NIR Grain Network Monthly Check Sample Results - HY2025-R02 – Wheat Flour (Spiked) – August 2025	5.6	4.92	-0.1



UK NIR Grain Network Monthly Check Sample Results - HY2025-R03 - Wheat Flour (Spiked) - August 2025	9.3	9.32	0.2
FAPAS Food Chemistry Proficiency Test - 04548 - Mycotoxins in Animal Feed (Cereal Based) - April - May 2025	46.8	46.0	-0.1
BIPEA 31B - Mycotoxins - Cereals - Rusk - 3 - 5431 - April 2025	7.88	7.28	-0.27
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx01] [Barley] - March 2025	4.50	4.54	0.02
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx02] [Wheat] - March 2025	0.10	<2.0	-
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx03] [Wheat] - March 2025	8.76	8.23	-0.18
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx01] [Wheat] - March 2025	10.52	9.68	-0.24
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx02] [Wheat] - March 2025	0.12	<2.0	-
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx03] [Wheat] - March 2025	0.12	<2.0	-
BIPEA 31E - Mycotoxins - Feed - Animal Feed - 15 - 3031 - January 2025	15.31	16.40	0.27



BIPEA 31B – Mycotoxins - Cereals – Maize Flour - 4 - 2231 – December 2024	8.94	7.92	-0.41
FAPAS Food Chemistry Proficiency Test - 04528 - Mycotoxins in Animal Feed (Cereal Based) - September – November 2024	191	181	-0.3
BIPEA 31B - Mycotoxins – Cereals – Wheat Flour –1- 6231 – October 2024	4.40	3.68	-0.55
FAPAS Food Chemistry Proficiency Test - 04519 – Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July – September 2024	15.3	16.4	0.3
FAPAS Food Chemistry Proficiency Test - 04520 - Multi - Mycotoxins in Maize Flour - July – September 2024	2.38	2.47	0.2
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Lentils - 2 - 5931 – May 2024	11.20	10.46	-0.24
FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April – May 2024	4.64	3.82	-0.8
BIPEA 31B – 7 - 188 - Mycotoxins – Cereals – Oats – 7 - 2931 – February 2024	6.65	5.96	-0.36
BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024	9.31	6.14	-1.23
FAPAS Food Chemistry Proficiency Test - 04498 – Mycotoxin Contamination in Animal Feed (Cereal Based) - November 2023 – January 2024	88.8	86.5	-0.1



BIPEA 31B - Mycotoxins - Cereals - Ground Corn - 6 - 187 - December 2023	Not Set	<1.5	-
FAPAS Food Chemistry Proficiency Test - 04482 - Multi - Mycotoxins in Maize Flour July - September 2023	2.52	2.88	0.7
FAPAS Food Chemistry Proficiency Test - 04481 - Mycotoxins in Pet Dog Food - July - August 2023	10.9	9.13	-0.8
BIPEA 31C - 20 - Mycotoxins - Baby Food - 21 - 3931 - June 2023	2.58	2.38	-0.26
BIPEA 31B - 185 Mycotoxin Contamination in Cereals - Rice - 06 - 4331 - June 2023	4.00	3.62	-0.32
BIPEA 31E - 3 - Mycotoxin Contamination in Animal feed - 13 - 3031 - January 2023	6.07	5.55	-0.29
BIPEA 31B - 183 Mycotoxin Contamination in Ground barley - 05 - 0531 - December 2022	9.08	7.97	-0.44
FAPAS Food Chemistry Proficiency Test - 17227 - Mycotoxins Ochratoxin A in Wine - August - September 2022	2.78	2.40	-0.6
FAPAS Food Chemistry Proficiency Test - 04449 - Mycotoxins in Dried Figs (Water/Fruit Slurry) - August - September 2022	10.4	9.45	-0.4
FAPAS Food Chemistry Proficiency Test - 04448 - Multi - Mycotoxins in Maize Flour - July - September 2022	2.72	3.38	1.1



FAPAS Food Chemistry Proficiency Test - 04439 - Multi - Mycotoxins in Maize Flour - April - May 2022	2.98	3.18	0.3
FAPAS Food Chemistry Proficiency Test - 04426 - Multi - Mycotoxins in Maize Flour - October - December 2021	3.31	3.40	0.1
FAPAS Food Chemistry Proficiency Test - 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021	3.79	3.60	-0.2
FAPAS Food Chemistry Proficiency Test - 04417 - Mycotoxins in Hazelnut (Water/Nut Slurry) - June - August 2021	7.35	7.93	0.4
FAPAS Food Chemistry Proficiency Test - 04411 - Multi - Mycotoxins in Maize Flour - April - June 2021	4.45	4.68	0.2
FAPAS Food Chemistry Proficiency Test - 04402 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2020 - January 2021	45.7	47.3	0.2
FAPAS Food Chemistry Proficiency Test - 04390 - Mycotoxins in Hazelnut (Water/Nut Slurry) - July - August 2020	7.96	6.87	-0.6
FAPAS Food Chemistry Proficiency Test - 04384 - Multi - Mycotoxins in Maize Flour - April - May 2020	4.82	6.19	1.3
FAPAS Food Chemistry Proficiency Test - 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	6.64	7.06	0.3