

Ring Test Results

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test 04378 Multi - Mycotoxins in Maize Flour February - March 2020	6.64	7.06	0.3
FAPAS Food Chemistry Proficiency Test 04384 Multi - Mycotoxins in Maize Flour April - May 2020	4.82	6.19	1.3
FAPAS Food Chemistry Proficiency Test 04390 Mycotoxins in Hazelnut (Water/Nut Slurry) July- August 2020	7.96	6.87	-0.6
FAPAS Food Chemistry Proficiency Test 04402 Mycotoxin Contamination in Animal Feed (Cereal Based) December 2020 - January 2021	45.7	47.3	0.2
FAPAS Food Chemistry Proficiency Test 04411 Multi - Mycotoxins in Maize Flour April - June 2021	4.45	4.68	0.2
FAPAS Food Chemistry Proficiency Test 04417 Mycotoxins in Hazelnut (Water/Nut Slurry) June - August 2021	7.35	7.93	0.4
FAPAS Food Chemistry Proficiency Test 04418 Multi - Mycotoxins in Maize Flour July - September 2021	3.79	3.60	-0.2



FAPAS Food Chemistry Proficiency Test 04426 Multi-Mycotoxins in Maize Flour October - December 2021	3.31	3.40	0.1
FAPAS Food Chemistry Proficiency Test 04439 Multi-Mycotoxins in Maize Flour April - May 2022	2.98	3.18	0.3
FAPAS Food Chemistry Proficiency Test 04448 Multi-Mycotoxins in Maize Flour July - September 2022	2.72	3.38	1.1
FAPAS Food Chemistry Proficiency Test 04449 Mycotoxins in Dried Figs (Water/Fruit Slurry) August - September 2022	10.4	9.45	-0.4
FAPAS Food Chemistry Proficiency Test 17227 Mycotoxins Ochratoxin A in Wine August - September 2022	2.78	2.40	-0.6