

### Ring Test Results

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test 04301 Multi Mycotoxins in Maize November-December 2016	3.03	3.28	0.4
BIPEA 31b 153 Mycotoxins - Cereals - Animal Feed - 09-3031 - February 2017	8.70	9.03	0.24
BIPEA 31b 154 Mycotoxins - Cereals - Bran - 02-3631 - April 2017	9.60	8.90	-0.26
FAPAS Food Chemistry Proficiency Test 04319 Multi Mycotoxins in Maize August - September 2017	4.73	5.40	0.6
FAPAS Food Chemistry Proficiency Test 04330 Multi Mycotoxins in Maize February - March 2018	2.75	2.75	0.0
FAPAS Food Chemistry Proficiency Test 04351 Mycotoxin Contamination in Animal Feed (cereal based) December 2018 - January 2019	30.2	28.2	-0.3
FAPAS Food Chemistry Proficiency Test 04354 Multi Mycotoxins in Maize February - March 2019	4.43	4.88	0.5
FAPAS Food Chemistry Proficiency Test 17192 Ochratoxin A in Dried Vine Fruit (Water/Fruit Slurry) April - May 2019	8.38	8.27	-0.1



## BIO-SHIELD

FAPAS Food Chemistry Proficiency Test 04359 Multi-Mycotoxins in Maize April - May 2019	2.69	2.74	0.1
FAPAS Food Chemistry Proficiency Test 04364 Mycotoxins in Paprika June - August 2019	22.3	20.2	-0.4
FAPAS Food Chemistry Proficiency Test 04366 Multi-Mycotoxins in Maize August - September 2019	1.99	2.04	0.1
FAPAS Food Chemistry Proficiency Test 04367 Mycotoxins in Dried Figs August- October 2019	4.75	5.23	0.5
FAPAS Food Chemistry Proficiency Test 04430 Mycotoxin Contamination in Animal Feed December 2021 - January 2022	74.50	66.85	-0.5
BIPEA 31E-2 Mycotoxin Contamination in Corn Silage 01-5831 May 2022	6.72	4.83	-0.98
BIPEA 31B-180 Mycotoxin Contamination in Rice 05-4331 June 2022	6.76	4.78	-1.02



Figure 1. Proficiency Test's z-Scores

