




**BIO-SHIELD T-2/HT-2 8**

Cat.Number B5648/B5696

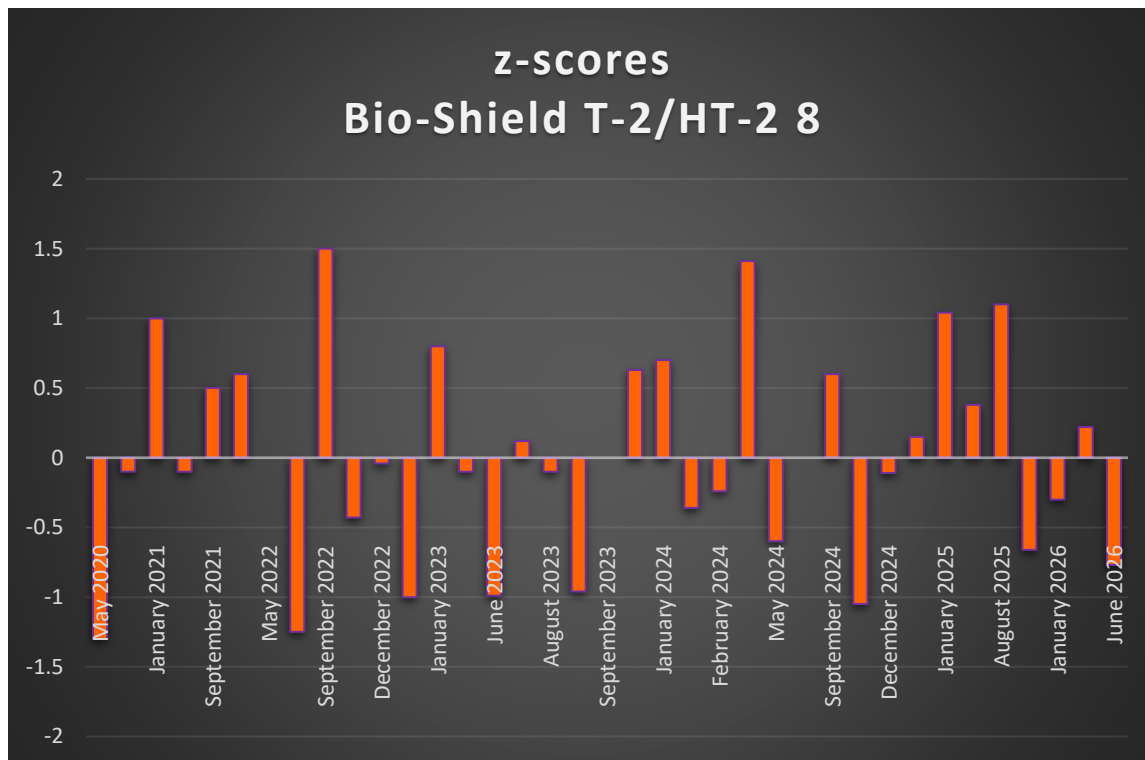

| Test                                                                                                           | Assigned value (µg/kg) | Result (µg/kg) | Z-score |
|----------------------------------------------------------------------------------------------------------------|------------------------|----------------|---------|
| BIPEA 31B - Mycotoxins – Cereals – Rice – 7 - 4331 – June 2026                                                 | 218                    | 175            | -0.77   |
| BIPEA 31E - Mycotoxins – Feed – Soyabean Meal – 18 - 6031 – May 2026                                           | 104                    | 110            | 0.22    |
| FAPAS Food Chemistry Proficiency Test - 04578 – Multi-mycotoxins in Durum Wheat - November 2025 – January 2026 | 118                    | 110            | -0.30   |
| BIPEA 31B - Mycotoxins – Cereals – Ground Corn – 19 - 0731 – December 2025                                     | 139                    | 115            | -0.66   |



|                                                                                                                             |      |       |       |
|-----------------------------------------------------------------------------------------------------------------------------|------|-------|-------|
| FAPAS Food Chemistry Proficiency Test - 04559 - Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July - August 2025 | 47.7 | 59.1  | 1.1   |
| BIPEA 31B - Mycotoxins - Cereals - Ground Rye - 4 - 5131 - June 2025                                                        | 35   | 39    | 0.38  |
| BIPEA 31E - Mycotoxins - Feed - Animal Feed - 15 - 3031 - January 2025                                                      | 48   | 62    | 1.04  |
| BIPEA 31C - Mycotoxins - Baby Food -24-3931 - December 2024                                                                 | 76   | 79    | 0.15  |
| BIPEA 31B - Mycotoxins - Cereals - Maize Flour - 4 - 2231 - December 2024                                                   | 102  | 99.0  | -0.11 |
| BIPEA 31B - Mycotoxins - Cereals - Wheat Flour -1- 6231 - October 2024                                                      | 115  | 83    | -1.05 |
| FAPAS Food Chemistry Proficiency Test - 04520 - Multi - Mycotoxins in Maize Flour - July - September 2024                   | 96.8 | 109   | 0.6   |
| BIPEA 31B - Mycotoxins - Cereals - Triticale - 3 - 1931 - June 2024                                                         | 79   | 70    | -0.42 |
| FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April - May 2024                        | 111  | 97.23 | -0.6  |
| BIPEA 31B - Cereals - Bran - 4 - 3631 - April 2024                                                                          | 102  | 140   | 1.41  |



|                                                                                                                  |      |       |       |
|------------------------------------------------------------------------------------------------------------------|------|-------|-------|
| BIPEA 31B – 7 - 188 - Mycotoxins – Cereals – Oats – 7 - 2931 – February 2024                                     | 223  | 209   | -0.24 |
| BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024                                              | 94   | 85    | -0.36 |
| FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 – January 2024 | 80.9 | 94.0  | 0.7   |
| BIPEA 31B - Mycotoxins – Cereals – Ground Corn - 6 - 187 – December 2023                                         | 210  | 244   | 0.63  |
| BIPEA 31B – 186 - Mycotoxin Contamination in Cereals – Common Wheat – 3 - 0631 - October 2023                    | 86.0 | 64.0  | -0.96 |
| FAPAS Food Chemistry Proficiency Test 04482 - Multi-Mycotoxins in Maize Flour - July - September 2023            | 104  | 103.2 | 0.0   |
| FAPAS Food Chemistry Proficiency Test 04481 - Mycotoxins in Pet Dog Food - July – August 2023                    | 40.2 | 39.1  | -0.1  |
| BIPEA 31c 20 Mycotoxins - Baby Food - 21 - 3931 – June 2023                                                      | 125  | 129   | 0.12  |
| BIPEA 31B - 185 Mycotoxin Contamination in Cereals – Rice - 06-4331 - June 2023                                  | 139  | 103   | -0.99 |
| FAPAS Food Chemistry Proficiency Test 04472 - Multi-Mycotoxins in Maize Flour - April – June 2023                | 169  | 166.7 | -0.1  |



|                                                                                                              |      |        |       |
|--------------------------------------------------------------------------------------------------------------|------|--------|-------|
| BIPEA 31B-183 Mycotoxin Contamination in Ground barley - 05-0531 - December 2022                             | 84.0 | 83.0   | -0.04 |
| BIPEA 31E-3 Mycotoxin Contamination in Animal feed - 13-3031 - January 2023                                  | 113  | 137    | 0.8   |
| FAPAS Food Chemistry Proficiency Test 22199 - Multi-Mycotoxins in Wheat Flour - November 2022-January 2023   | 93.7 | 72.54  | -1.0  |
| BIPEA 31B-182 Mycotoxin Contamination in Ground corn - 18-0731 - December 2022                               | 116  | 103    | -0.43 |
| FAPAS Food Chemistry Proficiency Test 04448 - Multi-Mycotoxins in Maize Flour - July - September 2022        | 92.6 | 123.42 | 1.5   |
| BIPEA 31B-180 Mycotoxin Contamination in Rice - 05-4331 - June 2022                                          | 99   | 66     | -1.25 |
| FAPAS Food Chemistry Proficiency Test 04439 - Multi-Mycotoxins in Maize Flour - April - May 2022             | 189  | 190.85 | 0.0   |
| FAPAS Food Chemistry Proficiency Test 22187 - Multi-Mycotoxins in Wheat Flour - November 2021 - January 2022 | 54   | 61.1   | 0.6   |
| FAPAS Food Chemistry Proficiency Test 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021      | 111  | 122.64 | 0.5   |
| FAPAS Food Chemistry Proficiency Test 04411 - Multi-Mycotoxins in Maize Flour - April - June 2021            | 174  | 169.47 | -0.1  |



|                                                                                                                                            |      |        |      |
|--------------------------------------------------------------------------------------------------------------------------------------------|------|--------|------|
| FAPAS Food Chemistry Proficiency Test<br>22177 - Multi - mycotoxins in Wheat Flour<br>- November 2020 - January 2021                       | 53.5 | 65.24  | 1.0  |
| FAPAS Food Chemistry Proficiency Test<br>22175 - Deoxynivalenol, Zearalenone, T-2<br>and HT-2 in Animal Feed - September -<br>October 2020 | 523  | 511.96 | -0.1 |
| FAPAS Food Chemistry Proficiency Test<br>04384 - Multi Mycotoxins in Maize Flour -<br>April - May 2020                                     | 330  | 247.37 | -1.3 |