


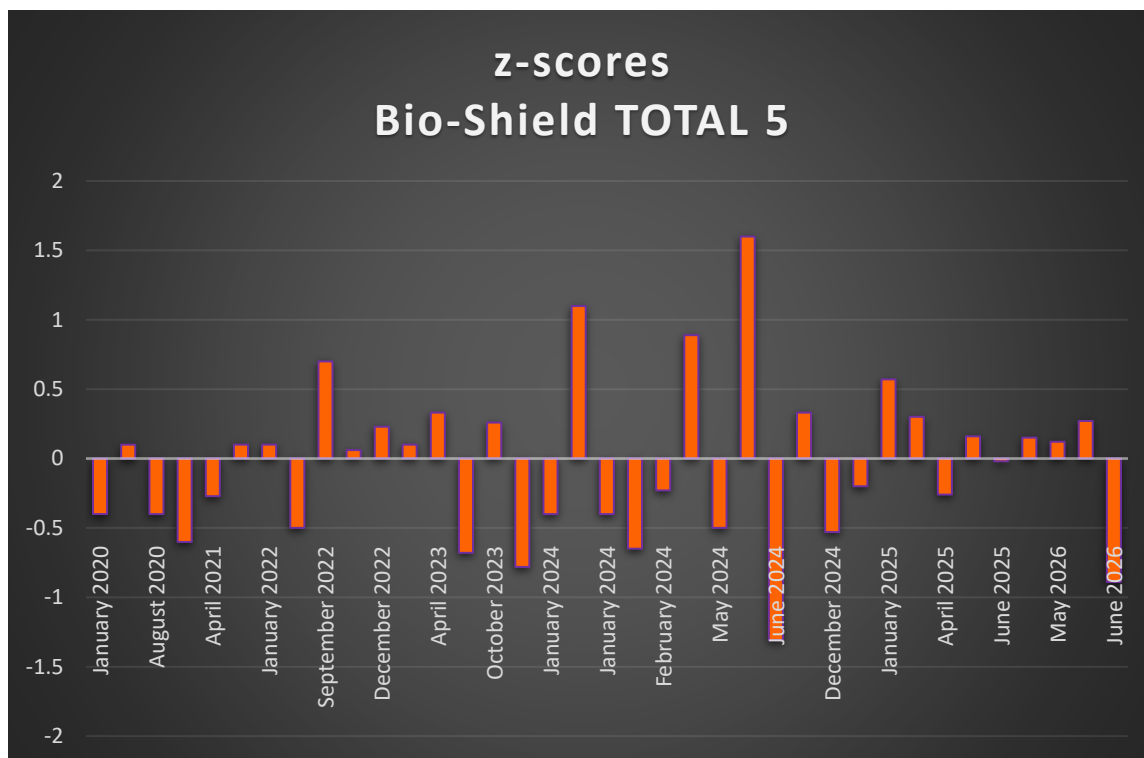

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31B - Mycotoxins – Cereals – Rice – 7 - 4331 – June 2026	15.70	12.00	-0.89
BIPEA 31E - Mycotoxins – Feed – Soyabean Meal – 18 - 6031 – May 2026	27.51	29.40	0.27
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Lentils – 03 - 5931 – May 2026	8.34	8.62	0.12
BIPEA 31B – Mycotoxins - Cereals – Pasta - 1 - 6431 – April 2026	3.95	4.13	0.15
BIPEA 31B – Mycotoxins - Cereals – Ground Rye - 4 - 5131 – June 2025	6.03	5.99	-0.02



BIPEA 31C – Mycotoxins – Baby Food - 25 - 3931 – June 2025	9.68	10.1	0.16
BIPEA 31B – Mycotoxins - Cereals – Rusk - 3 - 5431 – April 2025	11.49	10.68	-0.26
BIPEA 31B – Mycotoxins - Cereals – Rusk - 3 - 5431 – April 2025	13.62	11.7	-0.53
FAPAS Food Chemistry Proficiency Test - 04540 – Mycotoxins in Maize Flour - January – March 2025	4.93	5.23	0.3
BIPEA 31E – Mycotoxins - Feed – Animal Feed - 15 - 3031 – January 2025	33.07	37.90	0.57
BIPEA 31B – Mycotoxins - Cereals – Maize Flour - 4 - 2231 – December 2024	13.62	11.7	-0.53
FAPAS Food Chemistry Proficiency Test - 22231 - Multi - Mycotoxins in Wheat Flour - November 2024 – January 2025	8.63	8.30	-0.2
BIPEA 31B - Mycotoxins – Cereals – Wheat Flour -1- 6231 – October 2024	12.59	13.7	0.33
BIPEA 31B - Mycotoxins – Cereals – Triticale - 3 - 1931 – June 2024	13.14	8.51	-1.31
BIPEA 31E - Mycotoxins – Feed – Soyabean Meal - 16 - 6031 – May 2024	12.75	18.26	1.60
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Lentils - 2 - 5931 – May 2024	14.41	12.49	-0.50
BIPEA 31B - Cereals – Bran - 4 - 3631 – April 2024	15.76	19.50	0.89



BIPEA 31B – 7 - 188 - Mycotoxins – Cereals – Oats – 7 - 2931 – February 2024	7.02	6.55	-0.23
BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024	17.86	14.8	-0.65
FAPAS Food Chemistry Proficiency Test - 22215 - Multi - Mycotoxins in Wheat Flour - November 2023 – January 2024	5.90	5.44	-0.4
BIPEA 31B - Mycotoxins – Cereals – Ground Corn - 6 - 187 – December 2023	19.43	15.44	-0.78
FAPAS Food Chemistry Proficiency Test - 04497 - Aflatoxins in Peanut Butter (Water Slurry) - November 2023 – January 2024	8.91	11.07	1.1
FAPAS Food Chemistry Proficiency Test - 04496 - Aflatoxins in Rice – November 2023 – January 2024	9.77	9.02	-0.4
BIPEA 31B – 186 - Mycotoxin Contamination in Cereals – Common Wheat – 3 - 0631 - October 2023	5.06	5.45	0.26
BIPEA 31B – 185 - Mycotoxin Contamination in Cereals – Rice – 06 -4331 - June 2023	13.56	11.08	-0.68
BIPEA 31B – 184 - Mycotoxin Contamination in Cereals - 02-5431 - April 2023	8.13	8.89	0.33
FAPAS Food Chemistry Proficiency Test - 22199 – Multi - Mycotoxins in Wheat Flour - November 2022 - January 2023	8.88	9.11	0.1
BIPEA 31B - 183 - Mycotoxin Contamination in Ground Barley – 05 - 0531 - December 2022	19.32	20.50	0.23



BIPEA 31B - 182 Mycotoxin Contamination in Ground Corn - 18 - 0731 - December 2022	13.30	13.50	0.06
FAPAS Food Chemistry Proficiency Test - 04449 - Mycotoxins in Dried Figs (Water/Fruit Slurry) - August - September 2022	9.34	10.77	0.7
FAPAS Food Chemistry Proficiency Test - 04438 - Aflatoxins in Animal Feed (Cereal Based) - April - May 2022	19.10	16.78	-0.5
FAPAS Food Chemistry Proficiency Test - 22187 - Multi - Mycotoxins in Wheat Flour - November 2021 - January 2022	6.69	6.79	0.1
FAPAS Food Chemistry Proficiency Test - 04417 - Mycotoxins in Hazelnut (Water/Nut Slurry) - June - August 2021	10.2	10.42	0.1
QUALINOVA ROUND : AL 21008 - Aflatoxins in dried fruit - April 2021	9.64	9.07	-0.27
FAPAS Food Chemistry Proficiency Test - 04408 - Aflatoxins in Peanut (Water/Nut Slurry) - March - April 2021	12.1	10.55	-0.6
FAPAS Food Chemistry Proficiency Test - 04390 - Mycotoxins in Hazelnut (Water/Nut Slurry) - July - August 2020	9.67	8.79	-0.4
FAPAS Food Chemistry Proficiency Test - 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	6.87	6.99	0.1
FAPAS Food Chemistry Proficiency Test - 04374 - Aflatoxins in Rice - November 2019 - January 2020	9.31	8.4	-0.4