




Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test - 04551 – Mycotoxin Contamination in Infant Food – April – June 2025	0.36	<0.5	-
FAPAS Food Chemistry Proficiency Test - 04543 – Mycotoxins in Cereal Snack - February – April 2025	4.19	4.07	-0.1
FAPAS Food Chemistry Proficiency Test - 04546 - Aflatoxins in Hemp (CBD) Leaf - November 2024 – January 2025	10.6	9.44	-0.5
BIPEA 31C - Mycotoxins – Baby Food –24-3931 – December 2024	14.50	16.50	0.52

BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04528 - Mycotoxins in Animal Feed (Cereal Based) - September - November 2024		10.2	11.7	0.6
BIPEA 31C - Mycotoxins - Baby Food -23 - 3931 - June 2024		39.63	24.80	-1.46
BIPEA 31A - 188 - Mycotoxins - Dried Fruits, Spices and Other Products - Hazelnut Paste - 3 - 2031 - March 2024		19.51	20.60	0.21
BIPEA 31A - 186 Mycotoxins - Dried Fruits, Spices and Other Products - Ground Ginger - 02 - 1231 - November 2023		20.96	21.20	0.04
BIPEA 31A - Mycotoxins - Dried Fruits, Spices and Other Products - Peanut Paste - 6 - 185 - September 2023		8.69	9.57	0.36
BIPEA 31c 20 Mycotoxins - Baby Food - 21 - 3931 - June 2023		14.69	17.7	0.77
BIPEA 31A - 184 Mycotoxins Dried Fruits, Spices and Other Products - Pistachio Paste - 08-1431 May 2023		21.16	17.10	-0.73
QUALINOVA FOOD Mycotoxins in Feed - AL 23003 - April 2023	Sample 1	<1.0	0.989	-
	Sample 2	6.42	6.31	-0.07
BIPEA 31E-3 Mycotoxin Contamination in Animal feed 13 - 3031 January 2023		11.32	9.57	-0.57
BIPEA 31B-180 Mycotoxin Contamination in Rice 05 - 4331 June 2022		9.62	9.37	-0.09

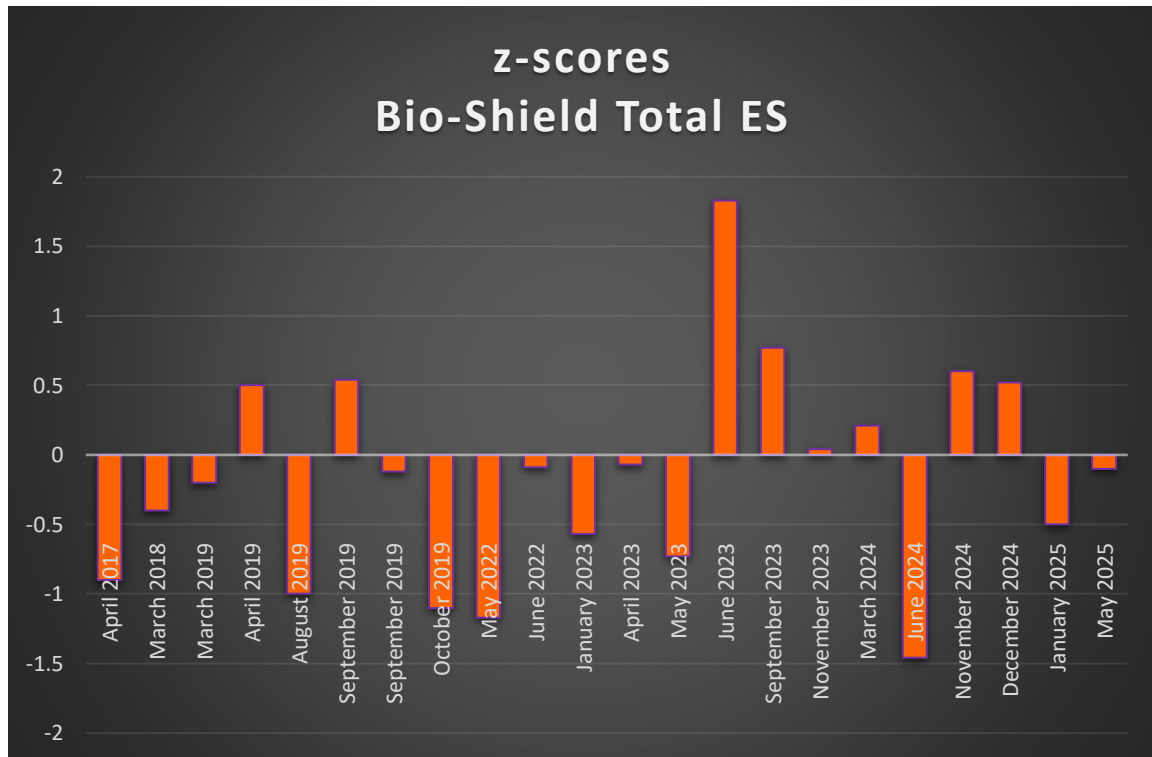

BIO-SHIELD

BIPEA 31E-2 Mycotoxin Contamination in Corn Silage 01 -5831 May 2022			6.48	4.29	-1.17
FAPAS Food Chemistry Proficiency Test 04367 Mycotoxins in Dried Figs August - October 2019			10.8	8.1	-1.1
DRRR 190663 AFLATOXIN 1	RVEP CORN	Sample 1	9.0	10.1	0.54
		Sample 2	17.9	16.6	-0.12
FAPAS Food Chemistry Proficiency Test 04364 Mycotoxins in Paprika June - August 2019			10.2	7.9	-1.0
FAPAS Food Chemistry Proficiency Test 04357 Aflatoxins in Peanut March - April 2019			12.40	13.86	0.5
FAPAS Food Chemistry Proficiency Test 04354 Multi Mycotoxins in Maize February - March 2019			4.95	4.78	-0.2
FAPAS Food Chemistry Proficiency Test 04330 Multi Mycotoxins in Maize February - March 2018			4.03	3.63	-0.4
BIPEA 31b 154 Mycotoxins - Cereals - Bran - 02-3631 - April 2017			28.0	21.5	-0.9



PROGNOSIS
BIOTECH