




Proficiency Test	Assigned value (mg/kg)	Result (mg/kg)	Z-score
FAPAS Egg in Cake Mix 27470AQC April – June 2026	5.81(allergenic protein)	5.23	-0.4
FAPAS Egg in Cake Mix 27470BQC April – June 2026	NOT DETECTED	<0.96	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1077 – June 2026	10.6	8.0	-0.90
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1177 – June 2026	18.5	19.5	0.59
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1277 – June 2026	NOT DETECTED	<0.9	-

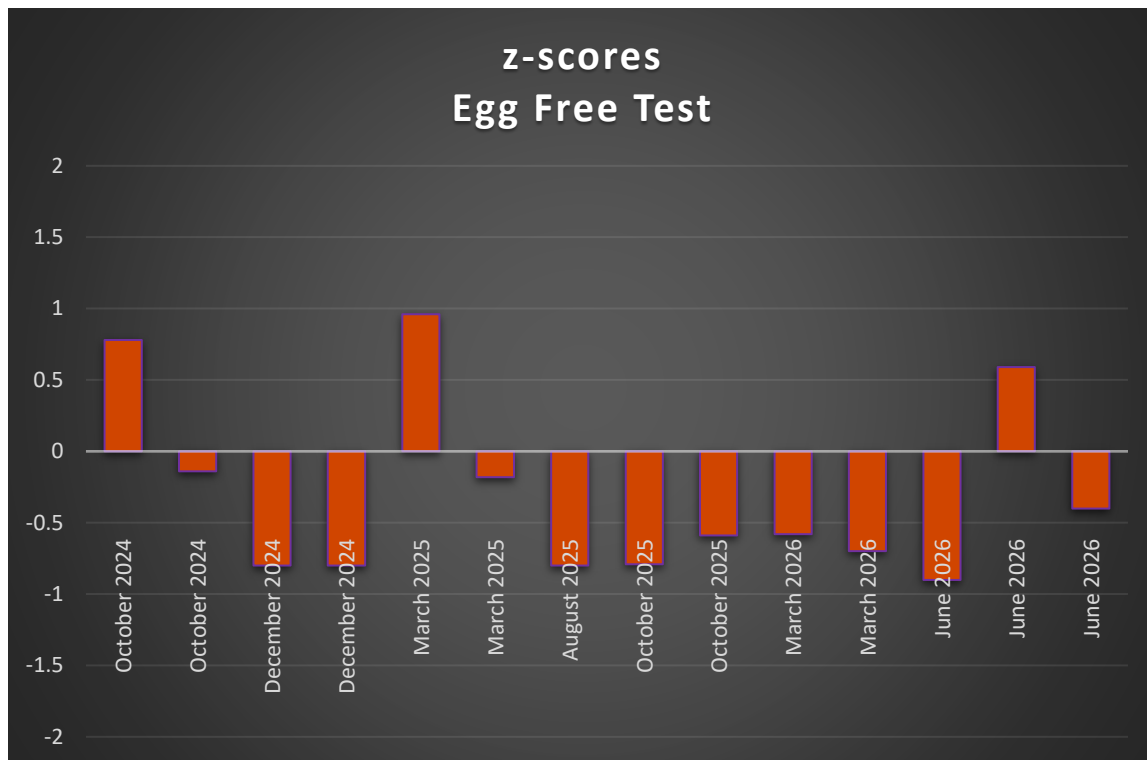

FAPAS Gluten, Milk & Egg in Cooked Biscuit 27459 January – March 2026	6.41	5.25	-0.7
BIPEA Food Allergens – EGG 77B – Pasta – 11 - 2577 - March 2026	42.5	33.0	-0.58
BIPEA Food Allergens – EGG 77B – Pasta – 10 - 2677 - March 2026	NOT DETECTED	<1.0	-
FAPAS Gluten, Milk & Egg in Cake Mix 27438B July – August 2025	18.8	15.07	-0.8
BIPEA Food Allergens – EGG 77B – Cake Mix – 11 - 0477 - October 2025	NOT DETECTED	<0.9	-
BIPEA Food Allergens – EGG 77B – Cake Mix – 11 - 0577 - October 2025	24.8	21.0	-0.59
BIPEA Food Allergens – EGG 77B – Cake Mix – 11 - 0677 - October 2025	45.4	39.5	-0.79
BIPEA Food Allergens – EGG 77B – Pasta – 10 - 2577 – March 2025	NOT DETECTED	<1.0	-
BIPEA Food Allergens – EGG 77B – Pasta – 9 - 2677 – March 2025	12.7	18.0	0.96
BIPEA Food Allergens – EGG 77B – Pasta – 9 - 2777 – March 2025	37.0	36.0	-0.18
FAPAS Gluten, Egg and Milk in Cake Mix 27400A July – September 2024	3.61	2.89	-0.8
FAPAS Gluten, Egg and Milk in Cake Mix 27400B July – September 2024	5.89	4.73	-0.8
BIPEA Food Allergens – EGG 77B – Cake Mix – 10 - 0677 - October 2024	23.5	22.4	-0.14
BIPEA Food Allergens – EGG 77B – Cake Mix – 10 - 0577 - October 2024	NOT DETECTED	<0.9	-
BIPEA Food Allergens – EGG 77B – Cake Mix – 10 - 0477 - October 2024	32.9	39.5	0.78



Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Allergens on Cotton Swabs 27440A August – October 2025	DETECTED	DETECTED	97.0
FAPAS Allergens on Cotton Swabs 27440B August – October 2025	NOT DETECTED	NOT DETECTED	97.0

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Egg Ei/Egg protein R6402	Range for $ z \leq 2$ according to R-Biopharm Ridascreen Fast Egg Ei/Egg protein R6402	Result (mg/kg)	Recovery (%)
FAPAS Gluten, Milk & Egg in Cake Mix 27438A July – August 2025	18.7	9.35 – 28.0	20.3	108.6
FAPAS Allergens in Sausage Meat (Raw) 27427A March – May 2025	16.8	8.42 – 25.3	22.04	131.2
FAPAS Allergens in Sausage Meat (Raw) 27427B March – May 2025	NOT DETECTED	-	<2.0	-
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27418 January – March 2025	5.33	2.67 – 8.00	2.80	52.5