




Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
FAPAS - Hazelnut in Chocolate – 27473A May – July 2026	NOT DETECTED	<2.0	-
FAPAS - Hazelnut in Chocolate – 27473B May – July 2026	3.32 (mg/kg of allergenic protein)	2.50	-1.0
BIPEA - Food Allergens - Nuts 77G - 21 - May 2026 - Cake - 14-2877	NOT DETECTED	<2.0	-
BIPEA - Food Allergens - Nuts 77G - 21 - May 2026 - Cake - 12-2977	59.0	70.0	0.75
BIPEA - Food Allergens - Nuts 77G - 21 - May 2026 - Cake - 16-3077	50.5	53.7	0.25

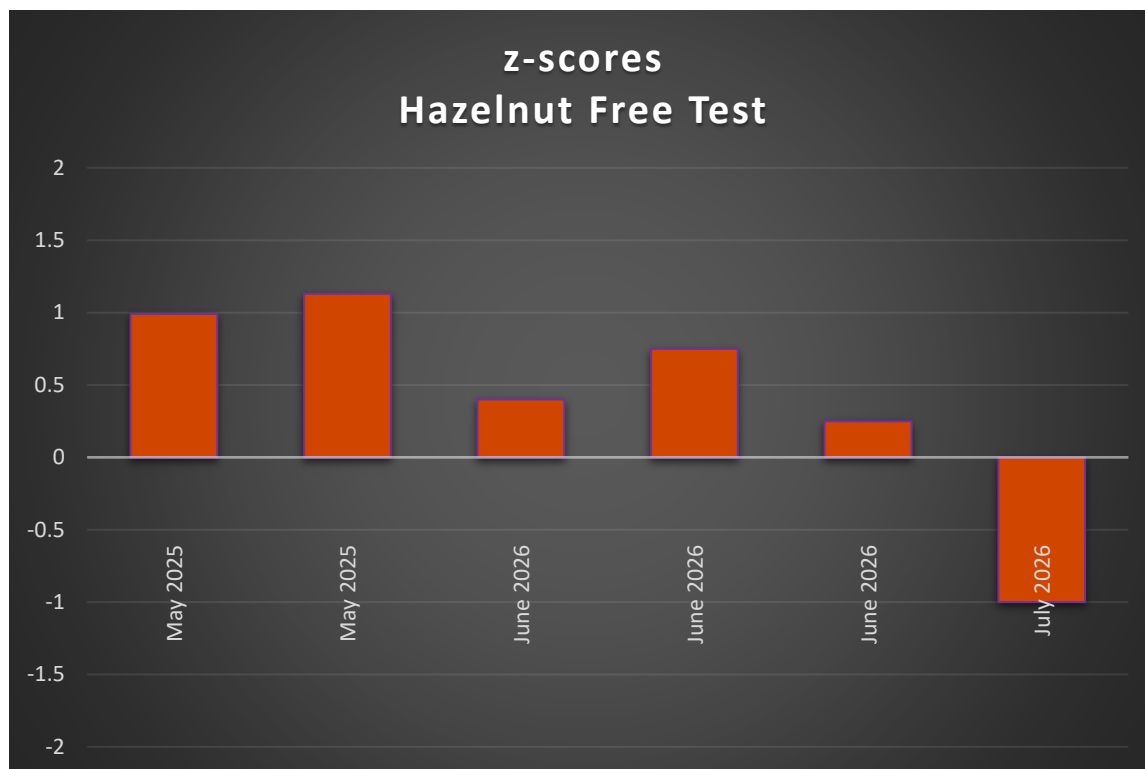

FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March – April 2026	3.69 (mg/kg of allergenic protein)	4.04	0.4
BIPEA - Food Allergens - Nuts 77G – 19 - May 2025 - Cake - 13-2877	532.9	705.5	0.99
BIPEA - Food Allergens - Nuts 77G – 19 - May 2025 - Cake - 11-2977	318.6	450.8	1.13
BIPEA - Food Allergens - Nuts 77G – 19 - May 2025 - Cake - 15-3077	NOT DETECTED	<2.0	-

Proficiency Test	Assigned value (mg/kg) according to ProGnosis Biotech Allergen-Shield Hazelnut A1648/A1696	Range for $ z \leq 2$ according to ProGnosis Biotech Allergen-Shield Hazelnut A1648/A1696	Result (mg/kg)	Recovery (%)
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March – April 2026	22.1	11.1 – 33.2	26.9	121.7

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS - Nuts In Chocolate - 27453 - November 2025 – January 2026	NOT DETECTED	<2.0	97.0
FAPAS - Hazelnut in Chocolate- 27433A - May – July 2025	DETECTED	DETECTED	100.0
FAPAS - Hazelnut in Chocolate- 27433B - May – July 2025	NOT DETECTED	<2.0	95.0



FAPAS - Hazelnut and Peanut in Cooked Biscuit -27425 - March - April 2025	DETECTED	DETECTED	95.0
FAPAS - Nuts In Chocolate - 27414 - November 2024 - January 2025	NOT DETECTED	NOT DETECTED	91.0
FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27389 - March - May 2024	DETECTED	DETECTED	96.0
FAPAS - Nuts In Chocolate - 27378 - November 2023 - January 2024	NOT DETECTED	<2.0	100.0