




Proficiency Test	Assigned value (mg/kg) of allergenic protein	Result (mg/kg)	Z-score
FAPAS Lupin in Flour 27465A March – May 2026	NOT DETECTED	<1.0	-
FAPAS Lupin in Flour 27465B March – May 2026	17.8	16.0	-0.4
FAPAS Lupin in Flour 27388A February – April 2024	19.6	17.5	-0.4
FAPAS Lupin in Flour 27388B February – April 2024	NOT DETECTED	<1.0	-

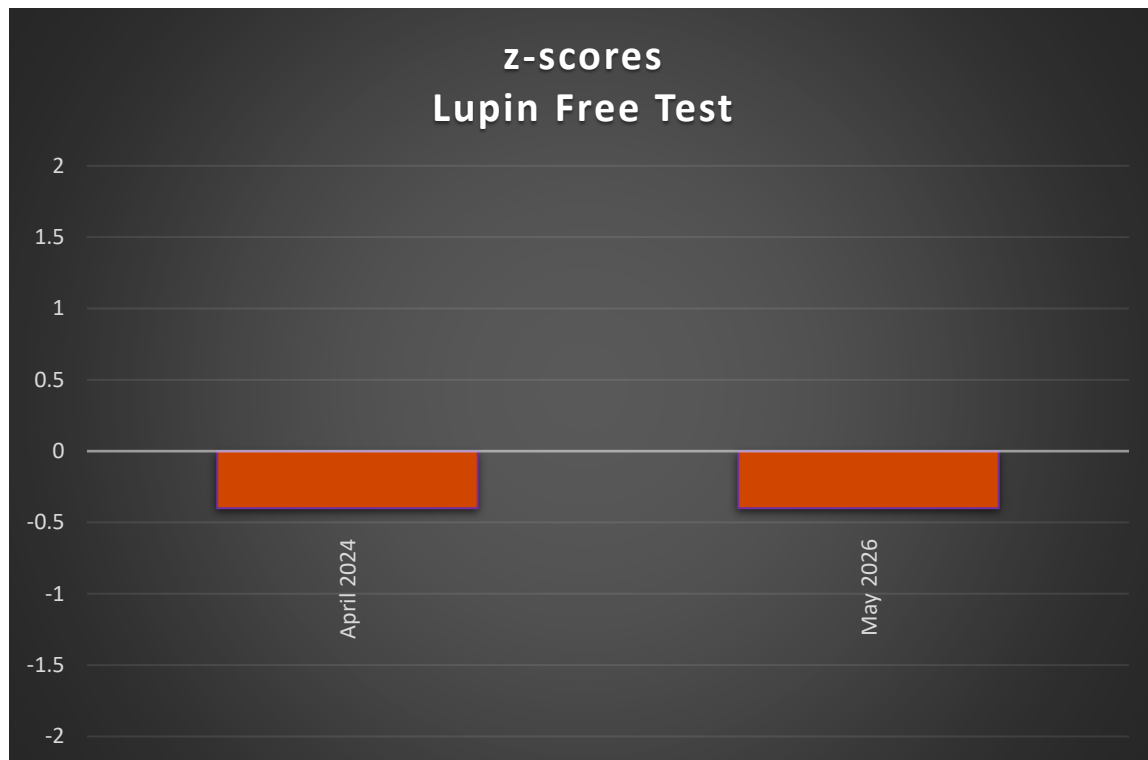

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm RIDASCREEN FAST Lupine R6102	Range for $ z \leq 2$ according to R-Biopharm RIDASCREEN FAST Lupine R6102	Result (mg/kg)	Recovery (%)
FAPAS Lupin in Flour 27424A March – April 2025	11.4	5.68 – 17.0	12.1	106.1
FAPAS Lupin in Flour 27424B March – April 2025	NOT DETECTED	-	<1.0	-
FAPAS Lupin in Flour 27388A February – April 2024	20.9	10.4 – 31.3	17.5	83.7
FAPAS Lupin in Flour 27388B February – April 2024	NOT DETECTED	-	<1.0	-