

| Proficiency Test  | Assigned value (mg/L)              | Result (mg/L) | Z-score |
|---|------------------------------------|---------------|---------|
| FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March - April 2026        | 5.26 (mg/kg of allergenic protein) | 7.25          | 1.5     |
| FAPAS Peanut as the Allergenic Protein in Chocolate 27437A - July - August 2025 | 3.76                               | 4.0           | 0.3     |
| FAPAS Peanut as the Allergenic Protein in Chocolate 27437B - July - August 2025 | NOT DETECTED                       | <0.5          | -       |

| Proficiency Test   | Assigned value (mg/kg) according to R-Biopharm Ridascreen Peanut R6811 | Range for $ z  \leq 2$ according to R-Biopharm Ridascreen Peanut R6811 | Result (mg/kg) | Recovery (%) |
|--|--|--|----------------|--------------|
| FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27466 March - April 2026               | 27.0   | 13.5 - 40.5  | 29.0           | 107.4        |
| FAPAS - Peanut as the Allergenic Protein in Chocolate - 27461A - February - March 2026 | NOT DETECTED   | -  | <2.0           | -            |
| FAPAS - Peanut as the Allergenic Protein in Chocolate - 27461B - February - March 2026 | 15.8   | 7.88 - 23.6  | 18.3           | 115.8        |
| FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27425 - March - April 2025             | 14.9   | 7.44 - 22.3  | 17.98          | 120.7        |
| FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27389 - March - May 2024               | NOT DETECTED   | -  | <2.0           | -            |

| Proficiency Test   | Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Peanut R6202 | Range for $ z  \leq 2$ according to R-Biopharm Ridascreen Fast Peanut R6202 | Result (mg/kg) | Recovery (%) |
|--|---|---|----------------|--------------|
| FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27425 - March - April 2025 | 34.5  | 17.3 - 51.8   | 22.6           | 65.5         |
| FAPAS - Hazelnut and Peanut in Cooked Biscuit - 27389 - March - May 2024   | 7.29  | 3.64 - 10.9   | 9.8            | 134.4        |

| Proficiency Test  | Assignment   | Result (mg/kg) | Satisfactory results / results agreeing with assignment (%) |
|---|--------------|----------------|---|
| FAPAS Peanut, Sesame & Soya on Swabs (Cotton) 27460A - January - March 2026 | NOT DETECTED | <2.0           | 100.0   |
| FAPAS Peanut, Sesame & Soya on Swabs (Cotton) 27460B - January - March 2026 | DETECTED     | DETECTED       | 90.0  |
| FAPAS - Nuts In Chocolate - 27453 - November 2025 - January 2026            | NOT DETECTED | <2.0           | 94.0  |
| FAPAS Allergens on Swab 27419A - May - July 2025                            | DETECTED     | DETECTED       | 100.0   |
| FAPAS Allergens on Swab 27419B - May - July 2025                            | NOT DETECTED | <0.5           | 92.0  |
| FAPAS - Peanut in Chocolate - 27420A - February - March 2025                | NOT DETECTED | <4.0           | 97.0  |
| FAPAS - Peanut in Chocolate - 27420B - February - March 2025                | DETECTED     | DETECTED       | 97.0  |
| FAPAS - Nuts In Chocolate - 27414 - November 2024 - January 2025            | NOT DETECTED | <4.0           | 97.0  |



|  |                 |          |      |
|--|-----------------|----------|------|
| FAPAS – Hazelnut and Peanut in Cooked Biscuit<br>-27389 March – May 2024 | DETECTED        | DETECTED | 94.0 |
| FAPAS Allergens on Swab 27383A - January -<br>March 2024                 | DETECTED        | DETECTED | 69.0 |
| FAPAS Allergens on Swab 27383B - January -<br>March 2024                 | NOT<br>DETECTED | <4.0     | 94.0 |