


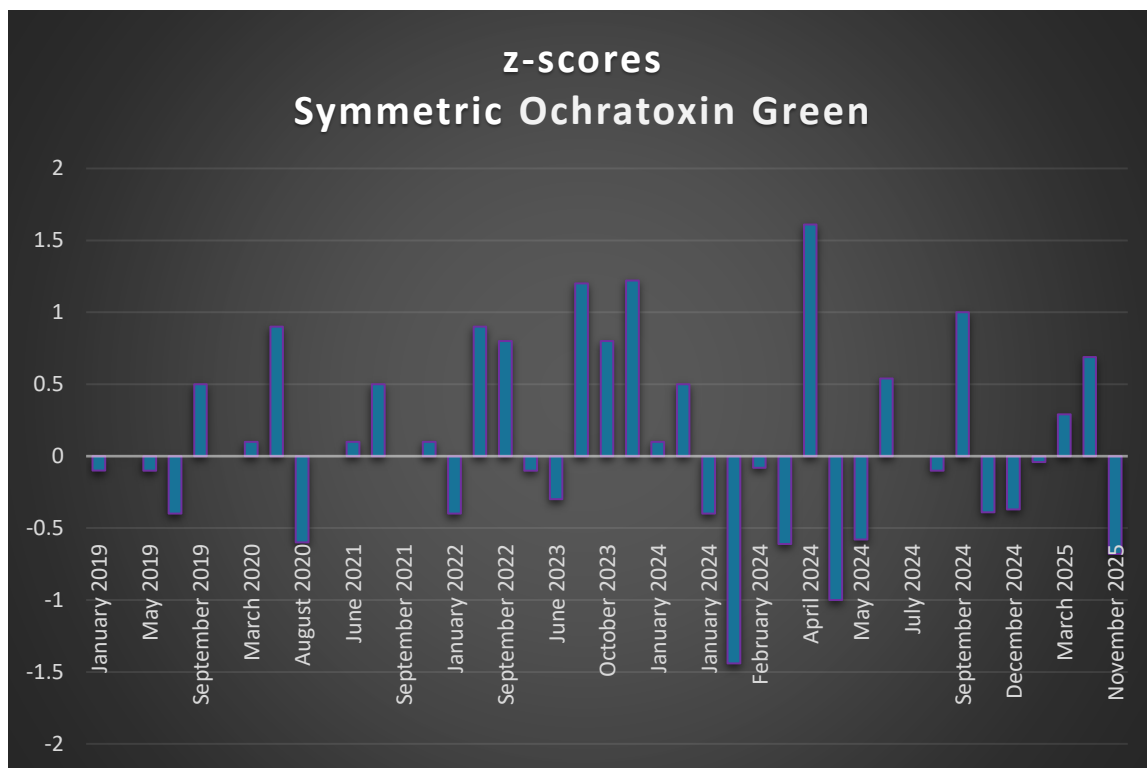

Test	Assinged value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Thyme – 01- 6331 – November 2025	8.05	6.50	-0.68
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx01] [Barley] – March 2025	4.50	4.41	-0.04
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R04-Tx02] [Wheat] – March 2025	0.10	<2.5	-
UK Grain Testing Network - Mycotoxins Ring Check Sample	8.76	9.64	0.29

[HY2024-R04-Tx03] [Wheat] - March 2025			
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx01] [Wheat] - March 2025	10.52	12.93	0.69
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx02] [Wheat] - March 2025	0.12	<2.5	-
UK Grain Testing Network - Mycotoxins Ring Check Sample [HY2024-R07-Tx03] [Wheat] - March 2025	0.12	<2.5	-
BIPEA 31B - Mycotoxins - Cereals - Maize Flour - 4 - 2231 - December 2024	8.94	8.02	-0.37
BIPEA 31B - Mycotoxins - Cereals - Wheat Flour -1- 6231 - October 2024	4.40	3.88	-0.39
FAPAS Food Chemistry Proficiency Test - 04519 - Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July - September 2024	15.3	18.5	1.0
FAPAS Food Chemistry Proficiency Test - 04520 - Multi - Mycotoxins in Maize Flour - July - September 2024	2.38	2.32	-0.1
BIPEA 31B - Mycotoxins - Cereals - Triticale - 3 - 1931 - June 2024	9.14	10.50	0.54

FAPAS Food Chemistry Proficiency Test - 04515 - Mycotoxins in Cocoa Powder – May – July 2024	4.56	4.57	0.0
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Lentils - 2 - 5931 – May 2024	11.20	9.44	-0.58
FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April – May 2024	4.64	3.59	-1.0
BIPEA 31B - Cereals – Bran - 4 - 3631 – April 2024	4.38	6.50	1.61
BIPEA 31A - 188 - Mycotoxins – Dried Fruits, Spices and Other Products – Hazelnut Paste – 3 - 2031 – March 2024	11.20	9.35	-0.61
BIPEA 31B – 7 - 188 - Mycotoxins – Cereals – Oats – 7 - 2931 – February 2024	6.65	6.50	-0.08
BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024	9.31	5.60	-1.44
BIPEA 31A – 5 - 187 - Mycotoxins – Dried Fruits, Spices and Other Products – Coffee – 5 - 1731 – January 2024	13.24	11.80	-0.40
FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 – January 2024	2.72	3.00	0.5
FAPAS Food Chemistry Proficiency Test - 04498 – Mycotoxin	88.8	90.0	0.1

Contamination in Animal Feed (Cereal Based) - November 2023 - January 2024			
BIPEA 31B - Mycotoxins - Cereals - Ground Corn - 6 - 187 - December 2023	Not Set	<1.0	-
BIPEA 31A - 186 Mycotoxins - Dried Fruits, Spices and Other Products - Ground Ginger - 02 - 1231 - November 2023	8.04	10.80	1.22
BIPEA 31B - 186 - Mycotoxin Contamination in Cereals - Common Wheat - 3 - 0631 - October 2023	5.25	6.50	0.8
FAPAS Food Chemistry Proficiency Test 04482 - Multi-Mycotoxins in Maize Flour - July - September 2023	2.52	3.20	1.2
FAPAS Food Chemistry Proficiency Test 04472 - Multi-Mycotoxins in Maize Flour - April - June 2023	4.12	3.89	-0.3
FAPAS Food Chemistry Proficiency Test 04449 - Mycotoxins in Dried Figs (Water/Fruit Slurry) - August - September 2022	10.4	10.12	-0.1
FAPAS Food Chemistry Proficiency Test 04448 - Multi-Mycotoxins in Maize Flour - July - September 2022	2.72	3.18	0.8
FAPAS Food Chemistry Proficiency Test 04439 - Multi-Mycotoxins in Maize Flour - April - May 2022	2.98	3.57	0.9

FAPAS Food Chemistry Proficiency Test 04430 - Mycotoxin Contamination in Animal Feed - December 2021 - January 2022	74.5	68.74	-0.4
FAPAS Food Chemistry Proficiency Test 04426 - Multi-Mycotoxins in Maize Flour - October - December 2021	3.31	3.42	0.1
FAPAS Food Chemistry Proficiency Test 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021	3.79	3.76	0.0
FAPAS Food Chemistry Proficiency Test 04417 - Mycotoxins in Hazelnut (Water/Nut Slurry) - June - August 2021	7.35	8.23	0.5
FAPAS Food Chemistry Proficiency Test 04411 - Multi-Mycotoxins in Maize Flour - April - June 2021	4.45	4.51	0.1
FAPAS Food Chemistry Proficiency Test 04402 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2020 - January 2021	45.7	45.18	0.0
FAPAS Food Chemistry Proficiency Test 04390 - Mycotoxins in Hazelnut (Water/Nut Slurry) - July - August 2020	7.96	6.85	-0.6
FAPAS Food Chemistry Proficiency Test 04384 - Multi - Mycotoxins in Maize Flour - April - May 2020	4.82	5.79	0.9
FAPAS Food Chemistry Proficiency Test 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	6.64	6.75	0.1

FAPAS Food Chemistry Proficiency Test 04367 - Mycotoxins in Dried Figs - August- October 2019	4.75	4.76	0.0
FAPAS Food Chemistry Proficiency Test 04366 - Multi-Mycotoxins in Maize - August - September 2019	1.99	2.21	0.5
FAPAS Food Chemistry Proficiency Test 04359 - Multi-Mycotoxins in Maize - April - May 2019	2.69	2.47	-0.4
FAPAS Food Chemistry Proficiency Test 17192 - Ochratoxin A in Dried Vine Fruit (Water/Fruit Slurry) - April – May 2019	8.38	8.17	-0.1
FAPAS Food Chemistry Proficiency Test 04354 - Multi Mycotoxins in Maize - February - March 2019	4.43	4.42	0.0
FAPAS Food Chemistry Proficiency Test 04351 - Mycotoxin Contamination in Animal Feed (cereal based) - December 2018 – January 2019	30.2	29.8	-0.1