




Proficiency Test	Assigned value (mg/kg)	Result (mg/kg)	Z-score
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1077 – June 2026	24.1	21.3	-0.44
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1177 – June 2026	NOT DETECTED	<3.0	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 21 - 1277 – June 2026	37.2	41.5	0.44
BIPEA Food Allergens - Milk 77C - Infant Formula – 23 - 0777 – April 2026	9.4	11.0	0.55
BIPEA Food Allergens - Milk 77C - Infant Formula – 23 - 0877 – April 2026	NOT DETECTED	<3.0	-
BIPEA Food Allergens - Milk 77C - Infant Formula – 22 - 0977 – April 2026	30.8	29.0	-0.23

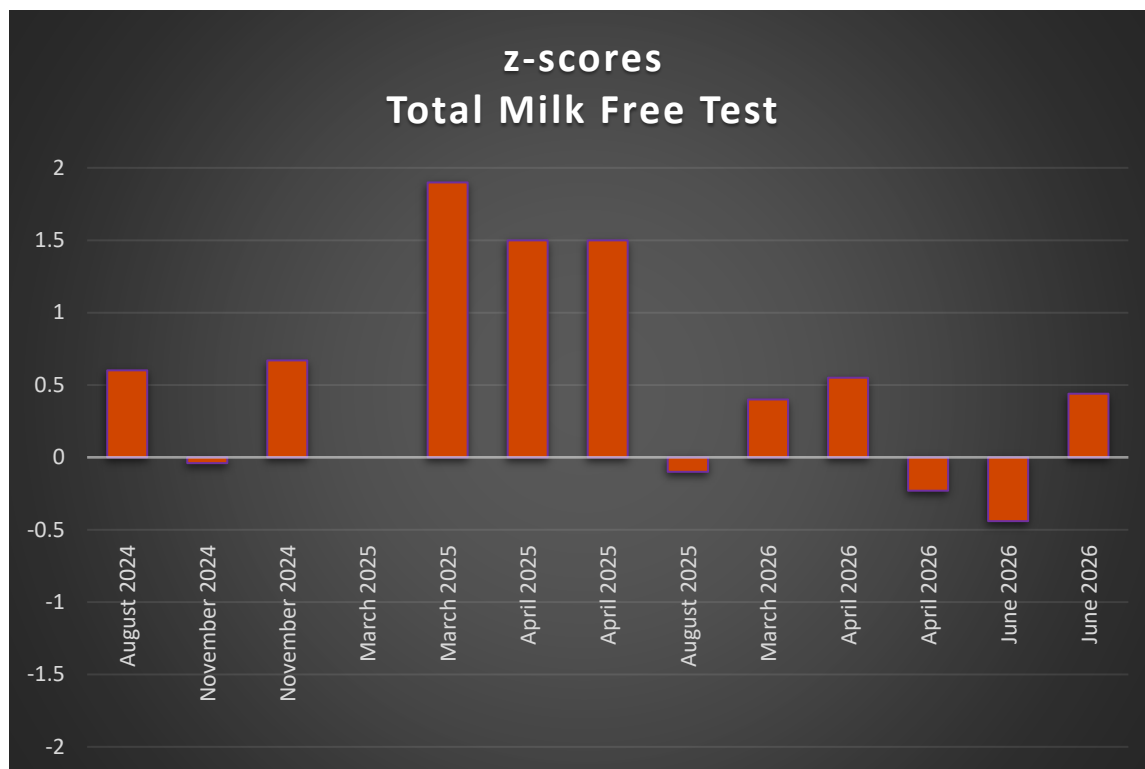

FAPAS Gluten, Milk & Egg in Cooked Biscuit 27459 January – March 2026	42.7	47.23	0.4
FAPAS Gluten, Milk & Egg in Cake Mix 27438B July – August 2025	62.0	60.0	-0.1
FAPAS Milk & Gluten in Plant Based Milk Alternative 27445A September – October 2025	NOT DETECTED	<4.0	-
FAPAS Milk & Gluten in Plant Based Milk Alternative 27445B September – October 2025	DETECTED	45.6	-
FAPAS Milk & Gluten in Infant Soya Formula 27422A February – March 2025	30.0	44.1	1.9
FAPAS Milk & Gluten in Infant Soya Formula 27422B February – March 2025	NOT DETECTED	<3.0	-
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 21 - 0777 – April 2025	NOT DETECTED	<3.0	-
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 21 - 0877 – April 2025	32.8	50.0	1.50
BIPEA Food Allergens - Milk 77C - 20 - Infant Formula – 20 - 0977 – April 2025	14.4	22.0	1.50
FAPAS Gluten, Milk & Egg in Cooked Biscuit 27418 January – March 2025	21.4	21.3	0.0
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 20 - 0777 – November 2024	NOT DETECTED	<3.0	-
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 20 - 0877 – November 2024	3.3	3.9	0.67
BIPEA Food Allergens - Milk 77C - 19 - Infant Formula – 19 - 0977 – November 2024	18.5	18.3	-0.04
FAPAS Milk in Infant Breakfast Cereal - 27398A - June – August 2024	NOT DETECTED	<2.5	-
FAPAS Milk in Infant Breakfast Cereal – 27398B - June – August 2024	33.8	38.4	0.60



Proficiency Test	Assigned value (mg/kg) according to Allergen-Shield Total Milk A1448/A1496	Range for $ z \leq 2$ according to Allergen-Shield Total Milk A1448/A1496	Result (mg/kg)	Recovery (%)
FAPAS Milk & Gluten in Infant Soya Formula 27463A March – May 2026	NOT DETECTED	-	<3.0	-
FAPAS Milk & Gluten in Infant Soya Formula 27463B March – May 2026	23.7	11.9 – 35.6	19.87	83.8

Proficiency Test	Assignment	Result (mg/kg)	Satisfactory results / results agreeing with assignment (%)
FAPAS Allergens on Cotton Swabs 27440A August – October 2025	NOT DETECTED	NOT DETECTED	95.0
FAPAS Allergens on Cotton Swabs 27440B August – October 2025	DETECTED	DETECTED	84.0
FAPAS Gluten, Egg and Milk in Cake Mix 27400A - July – September 2024	DETECTED	DETECTED	100%
FAPAS Gluten, Egg and Milk in Cake Mix 27400B - July – September 2024	DETECTED	DETECTED	100%

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Milk R4652	Range (mg/kg) for $ z \leq 2$ according to R-Biopharm Ridascreen Fast Milk R4652	Result (mg/kg)	Recovery (%)
FAPAS Gluten, Milk & Egg in Cake Mix 27438A July – August 2025	37.5	18.7 – 56.2	40.88	109.0
FAPAS Milk in Infant Breakfast Cereal 27435A June – August 2025	25.9	12.9 – 38.8	30.6	118.1



FAPAS Milk in Infant Breakfast Cereal 27435B June – August 2025	NOT DETECTED	-	<4.0	-
--	--------------	---	------	---